

HAPSITE® CDT Software and CDT IQ Laptop Software Installation Guide

Introduction

This document provides instructions for installing the HAPSITE CDT analytical module software and the CDT IQ laptop software.

These instructions assume the user has had HAPSITE CDT training and is familiar with the HAPSITE CDT analytical module and CDT IQ laptop software communications setup.

CDT IQ laptop software runs on Windows® 10 or 11 operating systems. CDT IQ software and software updates are available for download from the INFICON website under **Service & Support >> Download Software**.

NOTICE

If multiple HAPSITE CDT and laptops are being used, all must be updated to their respective software versions to maintain cross-system compatibility.

CAUTION

This software is designed only for HAPSITE CDT. Do not install this software on any other HAPSITE systems.

Installation of HAPSITE CDT Analytical Module Software


- 1 Confirm that the latest version of the HAPSITE CDT software has been downloaded from the website.
- 2 Open CDT IQ. Connect the laptop via wireless or Ethernet connection to the system. Confirm system is not preparing a method or in extended standby.




[ANALYZE] Heaters Not At Temperature

| Analyze: Cartridge_HT | | | | | NOT READY |
|-----------------------|-------------------------------------|-------------------------|-------------------------|------------------------|-----------|
| Target Chemicals | Detection Sensitivity Unknown | Collect Time 0:26 | Running Time 5:57 | Quant Library No | |
| | PREPARE ANALYZE | VIEW RESULTS | SELECT METHOD | EXTENDED STANDBY | |


- 3 In CDT IQ, right-click on the HAPSITE CDT sensor icon and select **Update HAPSITE Software**.

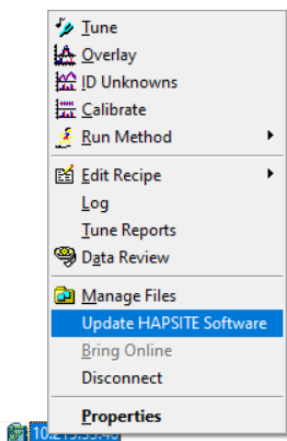


- Navigate to the location of the CDT SW download folder, select the file, and click **Open**.

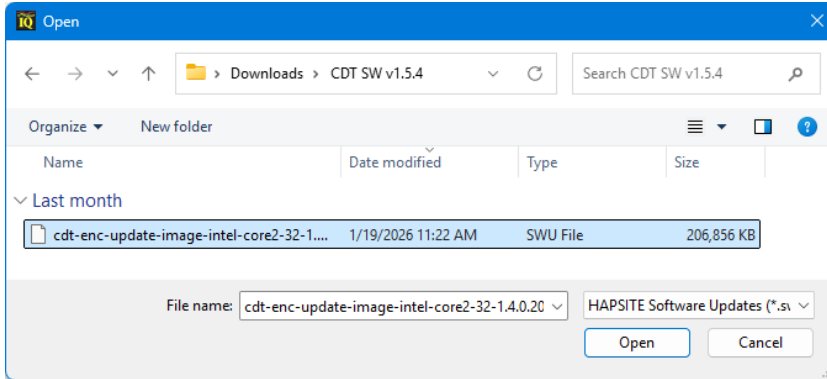


- A series of progress windows will indicate that the installation is in progress.

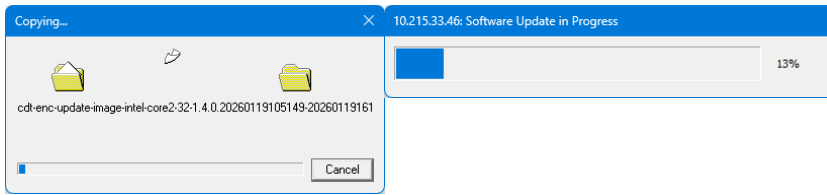

- A dialog box will be displayed upon completion. Click **Reboot** and allow system to reboot.

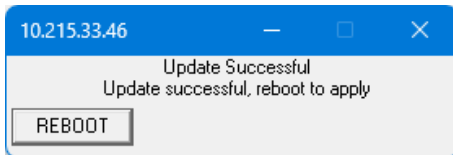


- Once the HAPSITE CDT system reboots, navigate to the System Settings screen on the front display. Confirm that the software has been updated.

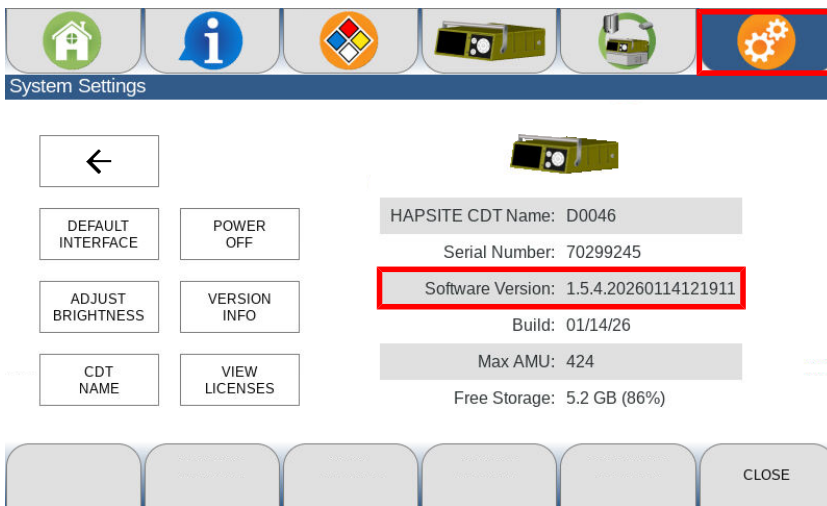


- Once confirmed, close out of CDT IQ. Proceed with updating the CDT IQ software.

Installation of CDT IQ Laptop Software


- 1 Navigate to the location of the CDT SW download folder. Open the CDT IQ setup file.

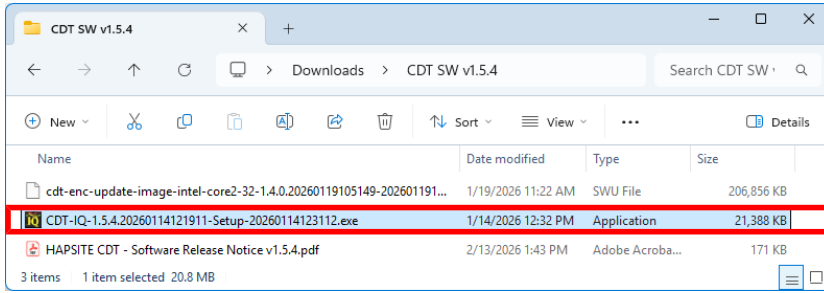


- 2 Follow the setup prompts. Click **Next**.

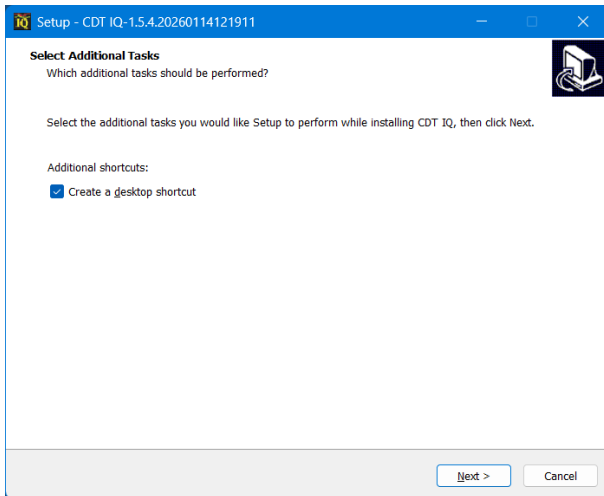


- 3 Click **Install**.

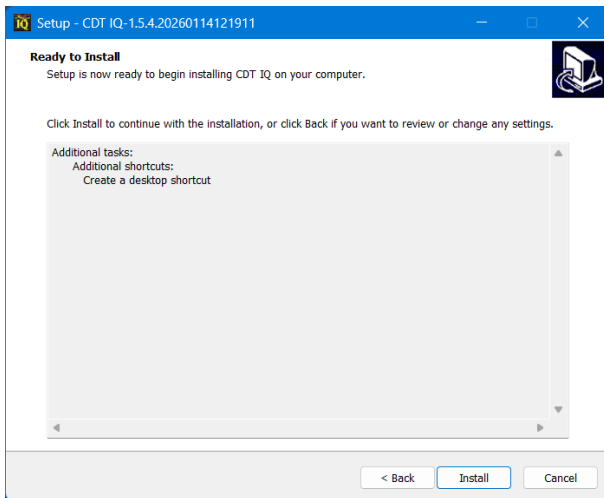

4 Allow software to complete installation.

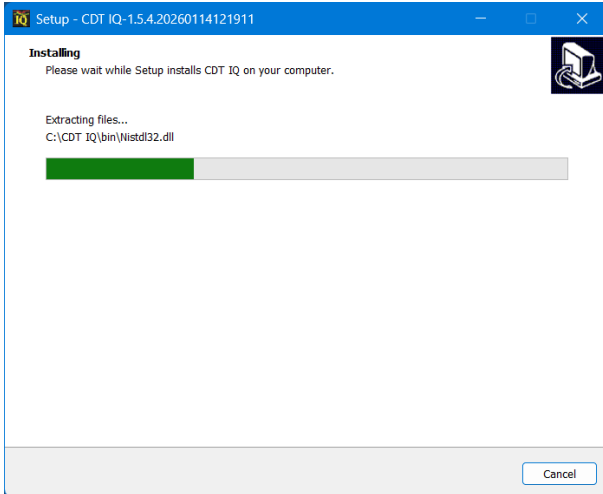

5 Click Finish.

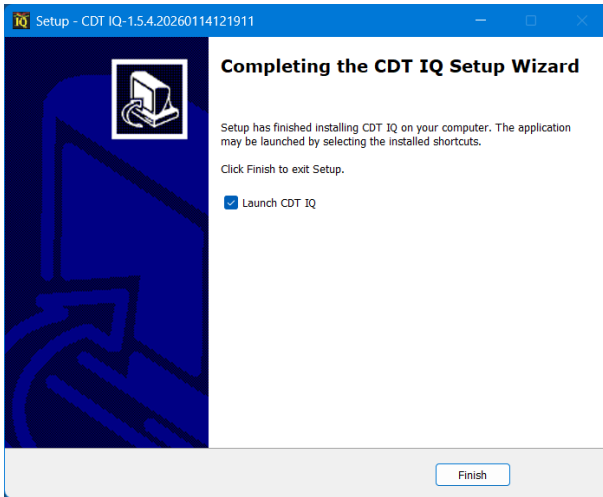

6 Open CDT IQ, click the Info icon, and confirm that the software has been updated.

