

HAPSITE® CDT

Software Release Notice v1.4.3

INFICON is pleased to release HAPSITE CDT Software v1.4.3.

Version 1.4.3 of the software is available for download from the INFICON website.

The latest release includes bug fixes and feature updates:

- A progress bar has been added during NEG activation.
- Lower MS dwell times are now supported, down to 200 μ s.
- A power button has been added to the touch display under the system settings screen.
- Data logging during methods has been added. Metadata (temperatures, pressures, etc) is saved to data file during methods. This data can be exported to .csv in CDT IQ.

Bug fixes

- Issues where the system would reboot instead of shutting down are addressed in this bug fix.
- Onboard AMDIS is now using the correct scan direction.
- Methods will now abort when accessories are uninstalled during runs. Additionally, to avoid the following popup error, please touch close after analysis is finished, prior to detaching an accessory.

