



3000 Micro GC

Gas Analyzer

PN 074-558-P1A

DETECT TO PROTECT™



R E F E R E N C E G U I D E

3000 Micro GC

Gas Analyzer

PN 074-558-P1A



www.inficon.com reachus@inficon.com

©2013 INFICON

Trademarks

The trademarks of the products mentioned in this manual are held by the companies that produce them.

3000 Micro GC® and Opening The Field To New Ideas® are registered trademarks of INFICON.

Windows®, and Microsoft® are registered trademarks of Microsoft Corporation.

Swagelok® is a registered trademark of Swagelok. All other brand and product names are trademarks or registered trademarks of their respective companies.

Disclaimer

The information contained in this manual is believed to be accurate and reliable. However, INFICON assumes no responsibility for its use and shall not be liable for any special, incidental, or consequential damages related to the use of this product.

Due to our continuing program of product improvements, specifications are subject to change without notice.

Copyright

©2013 All rights reserved.

Reproduction or adaptation of any part of this document without permission is unlawful.

Purpose of this Reference Guide

3000 Micro GC Guides are not intended to replace the 3000 Micro GC Operating Manual (PN 074-519-P1) or 3000 Micro GC training, but are intended for use as supporting materials for those who have been trained in an INFICON® sanctioned training course.

Definition of Note, Hint, Danger, Warning and Caution Paragraphs

NOTE: These messages contain pertinent information that is useful for achieving maximum 3000 Micro GC efficiency.

HINT: These messages provide insight into product usage.



DANGER

Failure to heed these messages may result in serious personal injury or death!



WARNING

Failure to heed these messages could result in personal injury.



WARNING - Risk Of Electric Shock

Dangerous voltages are present. Failure to heed these messages could result in personal injury.



CAUTION

Failure to heed these messages could result in damage to 3000 Micro GC.

How To Contact Customer Support

Worldwide support information regarding:

- ♦ Technical Support, to contact an applications engineer with questions regarding INFICON products and applications,
or
- ♦ Sales and Customer Service, to contact an INFICON Sales office,
or
- ♦ Repair Service, to contact an INFICON Service Center, is available at www.inficon.com.

If experiencing an instrument problem, please have the following information readily available:

- ♦ the instrument serial number,
- ♦ a description of the problem,
- ♦ an explanation of any corrective action that may have already been attempted,
- ♦ and the exact wording of any error messages observed.

To contact Customer Support, see Support at www.inficon.com.

This page is intentionally blank.

Table Of Contents

Cover Page	
Title Page	
Trademarks	
Disclaimer	
Copyright	
Purpose of this Reference Guide	
Definition of Note, Hint, Danger, Warning and Caution Paragraphs	
How To Contact Customer Support	
3000 Micro GC Guide 1	
Configuring a PC for a Direct Connection to	
3000 Micro GC via the Crossover Cable	1-1
3000 Micro GC Guide 2	
Connecting to 3000 Micro GC	
via the Static IP Address	2-1
3000 Micro GC Guide 3	
Changing the Instrument IP Address	3-1
3000 Micro GC Guide 4	
Changing the Carrier Gas Configuration	4-1
3000 Micro GC Guide 5	
Installing EZ IQ	5-1
3000 Micro GC Guide 6	
Configuring a 3000 Micro GC with EZ IQ	6-1
3000 Micro GC Guide 7	
3000 Micro GC Method Setup	7-1
3000 Micro GC Guide 8	
Running a Sample	8-1
3000 Micro GC Guide 9	
Integrating Chromatograms	9-1

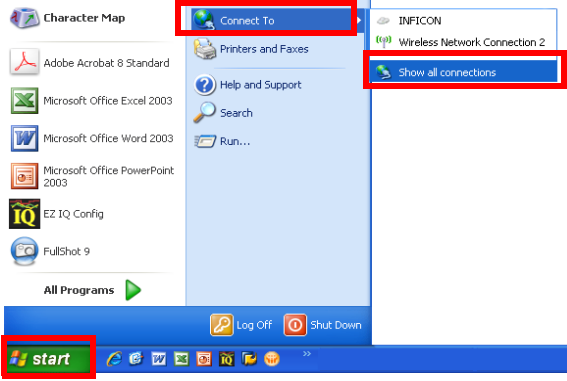
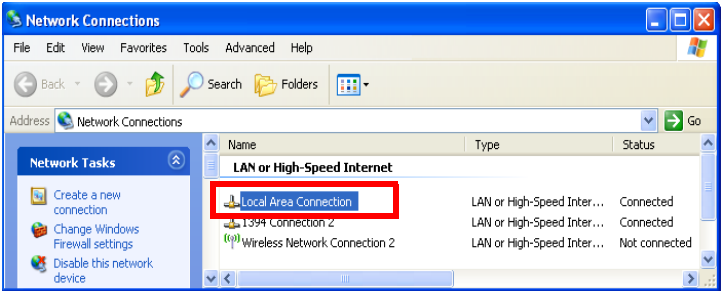
3000 Micro GC Guide 10 Setting Annotations	10-1
3000 Micro GC Guide 11 Qualifying Peaks and Groups	11-1
3000 Micro GC Guide 12 Adjusting Retention Time Windows	12-1
3000 Micro GC Guide 13 Conducting a Single Level Calibration	13-1
3000 Micro GC Guide 14 Conducting a Multiple Level Calibration	14-1
3000 Micro GC Guide 15 Calibration with Average Replicates	15-1
3000 Micro GC Guide 16 Customizing a Method Report	16-1
3000 Micro GC Guide 17 Customizing an ESTD Report	17-1
3000 Micro GC Guide 18 Changing a Standard Module	18-1
3000 Micro GC Guide 19 Changing a Performance Enhanced Module	19-1
3000 Micro GC Guide 20 Replacing a Standard Module with a Performance Enhanced Module	20-1
3000 Micro GC Guide 21 Testing Instrument Flow Rates	21-1
3000 Micro GC Guide 22 Installing the Digital I/O Module	22-1
3000 Micro GC Guide 23 Replacing the Battery Cable Fuse on a Portable Instrument	23-1

3000 Micro GC Guide 1

Configuring a PC for a Direct Connection to

3000 Micro GC via the Crossover Cable

Table 1-1 Configuring a PC for a direct connection via the crossover cable

Step	Description
1	Connect one end of the crossover cable (PN 5183-4649 provided in the ship kit) to a TCP/IP ethernet port on the PC. Connect the other end to the ethernet port on the back of 3000 Micro GC.
2	Turn the 3000 Micro GC power switch, located on its front, to the on position. The switch will illuminate green to indicate that 3000 Micro GC is on. Wait three minutes for the instrument to load.
3	<p>Click (Windows) Start >> Connect to >> Show All Connections.</p> <p>NOTE: Depending on the version of Windows, this navigation path may vary.</p>  A screenshot of the Windows XP Start menu. The 'Start' button is highlighted with a red box. The 'Connect To' menu item is highlighted with a red box. The 'Show all connections' option in the 'Connect To' submenu is also highlighted with a red box.
4	<p>On the Network Connections window, double-click Local Area Connection.</p>  A screenshot of the Windows 'Network Connections' window. The 'Local Area Connection' link is highlighted with a red box. The window shows a list of network connections with columns for Name, Type, and Status. The connections listed are 'LAN or High-Speed Internet', 'Local Area Connection', '1394 Connection 2', and 'Wireless Network Connection 2'.

PN 074-558-P1A

Table 1-1 Configuring a PC for a direct connection via the crossover cable (continued)

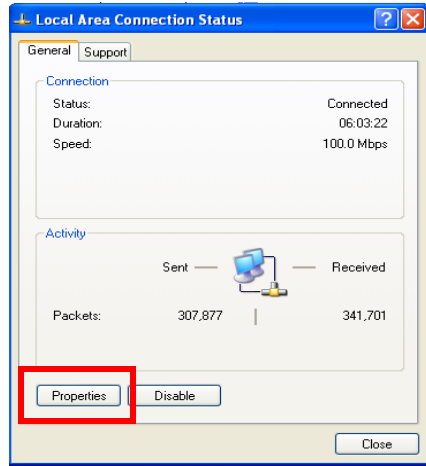
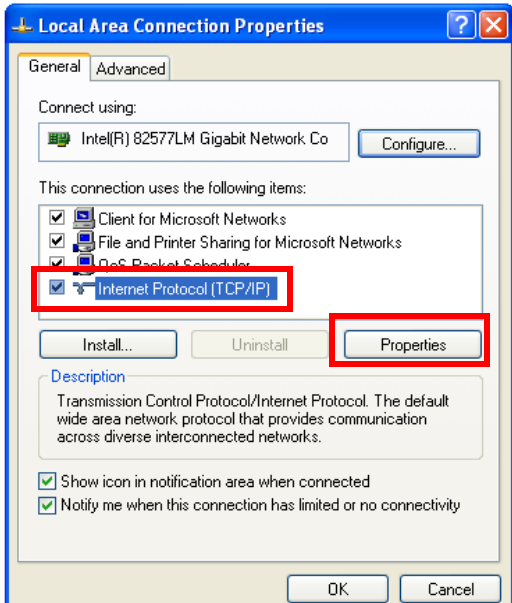
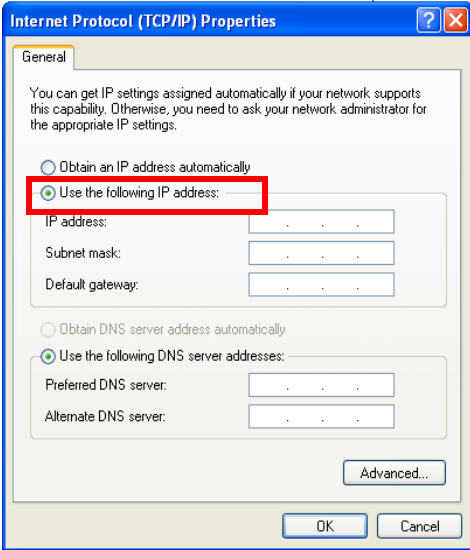
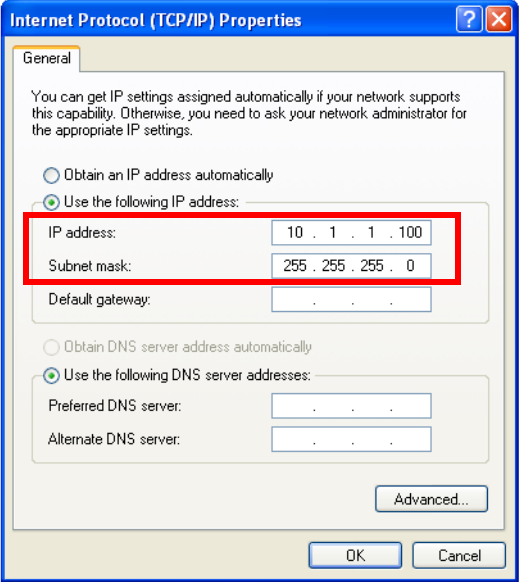

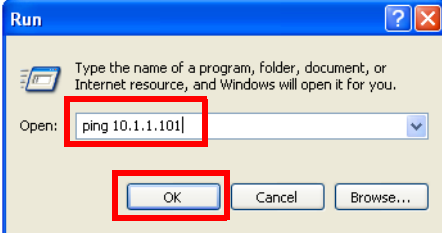
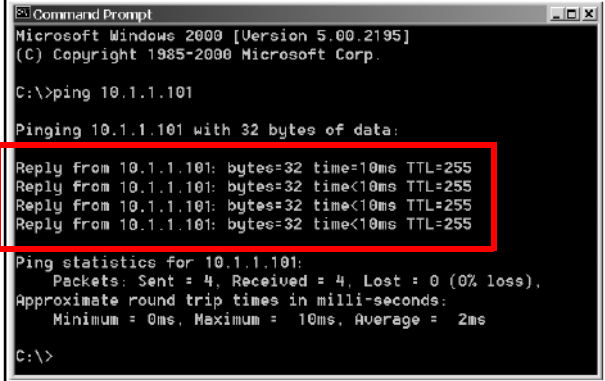
Step	Description
5	<p>In the Local Area Connection Status window, click Properties.</p> 
6	<p>In the Local Area Connection Properties window in The connection uses the following items: area, select Internet Protocol (TCP/IP), then click Properties.</p> 

Table 1-1 Configuring a PC for a direct connection via the crossover cable (continued)

Step	Description
7	<p>In the Internet Protocol (TCP/IP) Properties window, select Use the following IP address.</p> 
8	<p>Type the PC IP address and Subnet mask as follows:</p> <p>IP address: 10.1.1.100</p> <p>Subnet mask: 255.255.255.0</p> <p>Default gateway: *</p> <p>* The Default gateway can remain blank for a direct connection via the crossover cable.</p> 
9	Click OK to exit the Internet Protocol (TCP/IP) Properties window.
10	Click OK to exit the Local Area Connection Properties window.
11	Click Close to exit the Local Area Connection Status window.

PN 074-558-P1A

Table 1-1 Configuring a PC for a direct connection via the crossover cable (continued)

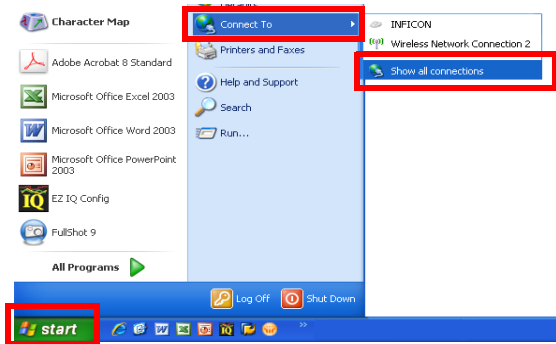
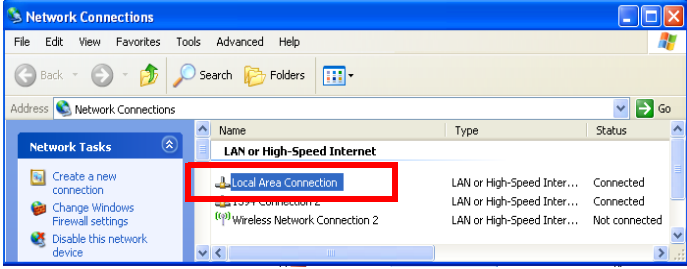
Step	Description
12	Click (Windows) Start >> Run . 
13	In the Run window, in the Open: box, type ping 10.1.1.101 . Click OK . 
14	Verify from the Command Prompt that the PC is receiving data packets from the instrument. NOTE: If the command prompt reads Request Timed Out or Detection Limit Unreachable , repeat steps 3 to 7 to verify IP settings. Verify that 3000 Micro GC is turned on and connected to the PC via the crossover cable. If there is still no communication contact INFICON. 

3000 Micro GC Guide 2

Connecting to 3000 Micro GC

via the Static IP Address

Table 2-1 Connecting to the instrument via the static IP address

Step	Description
1	<p>3000 Micro GC has both a static IP address and a dynamic IP address. The factory default dynamic IP address is set at 10.1.1.101. In the event that the IP address is modified or lost, it is possible to access the instrument through its static IP address: 192.168.1.99.</p> <p>NOTE: It is necessary to access the IP address via a direct crossover cable, in order to avoid potential routing issue in a local area network (LAN).</p>
2	<p>Connect one end of the crossover cable (PN 5183-4649 provided in the ship kit) to a TCP/IP ethernet port on the PC. Connect the other end to the ethernet port on the back of 3000 Micro GC.</p>
3	<p>The 3000 Micro GC power switch is located on its front panel. Turn the switch to the on position. The switch will illuminate green to indicate that 3000 Micro GC is on. Wait three minutes for the instrument to load.</p>
4	<p>Click (Windows) Start >> Connect to >> Show All Connections.</p> <p>NOTE: Navigation path may vary, depending on what version of Windows is running.</p> 
5	<p>On the Network Connections window, double-click Local Area Connection.</p> 

PN 074-558-P1A

Table 2-1 Connecting to the instrument via the static IP address (continued)

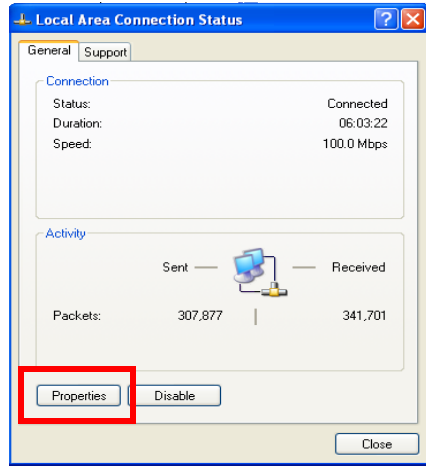
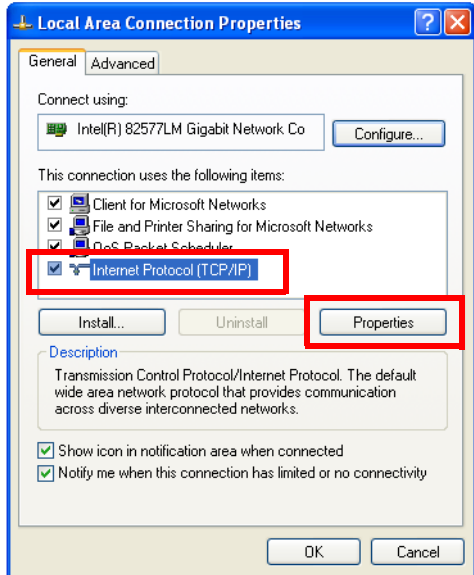
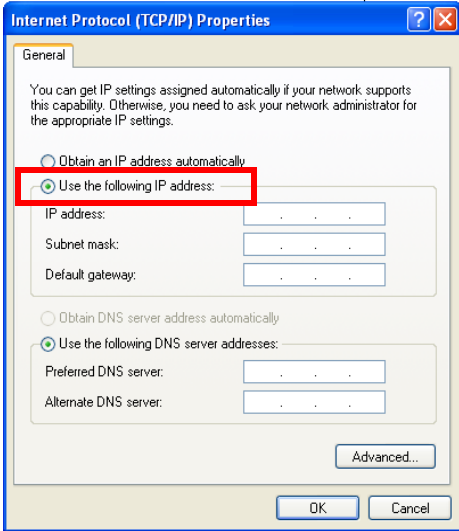
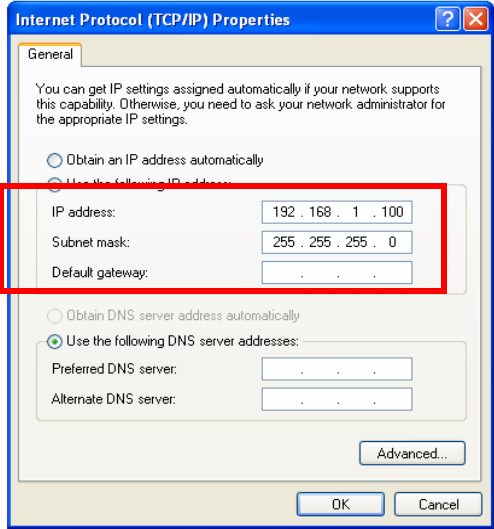

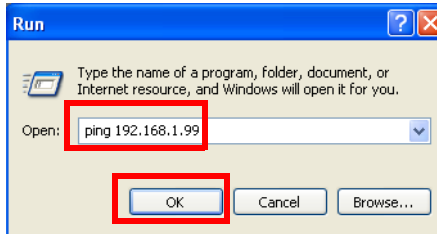
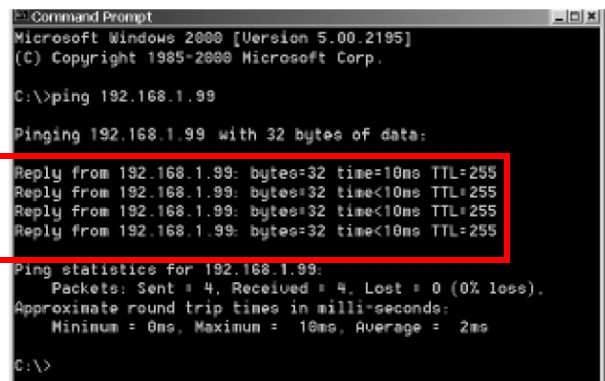
Step	Description
6	<p>In the Local Area Connection Status window, click Properties.</p> 
7	<p>The Local Area Connection Properties window will display. Select Internet Protocol (TCP/IP) in The connection uses the following items: area, then click Properties.</p> 

Table 2-1 Connecting to the instrument via the static IP address (continued)

Step	Description
8	<div>In the Internet Protocol (TCP/IP) Properties window, select Use the following IP address.</div> <div></div>
9	<div>Type the PC IP Address and Subnet Mask as follows:</div> <div>IP Address: 192.168.1.100</div> <div>Subnet Mask: 255.255.255.0</div> <div>Default gateway: *</div> <div>* The Default gateway can remain blank for a direct connection via the crossover cable.</div> <div></div>
10	Click OK to exit the Internet Protocol (TCP/IP) Properties window.
11	Click OK to exit the Local Area Connection Properties window.
12	Click Close to exit the Local Area Connection Status window.

PN 074-558-P1A

Table 2-1 Connecting to the instrument via the static IP address (continued)

Step	Description
13	Click (Windows) Start >> Run . 
14	In the Run window, in the Open: box, type ping 192.168.1.99 . Click OK . 
15	Verify from the Command Prompt that the PC is receiving data packets from the instrument. NOTE: If the command prompt reads Request Timed Out or Detection Limit Unreachable , repeat steps 3 to 7 to verify IP settings. Verify that 3000 Micro GC is on connected to the PC via the crossover cable. If there is still no communication contact INFICON. 
16	Open a web browser, such as Internet Explorer.
17	In the address bar, type 192.168.1.99 and press Enter .
18	The instrument embedded web page will display. Select the IP Configuration tab. Review the IP address.

3000 Micro GC Guide 3

Changing the Instrument IP Address

Table 3-1 Changing the Instrument IP Address

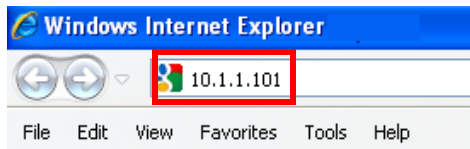
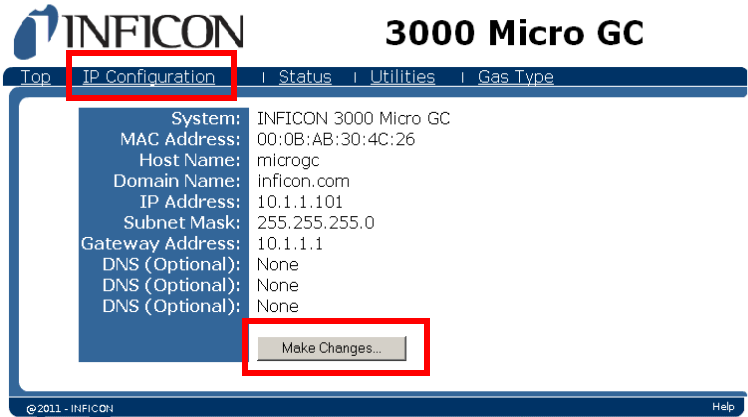
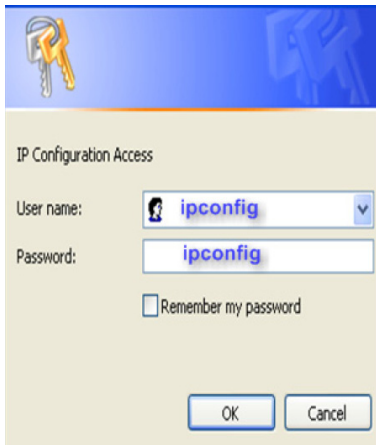
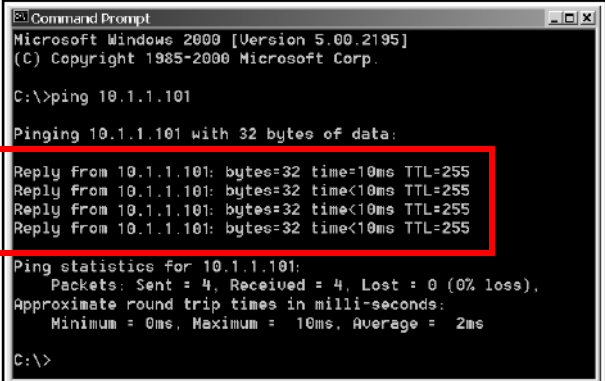
Step	Description
1	<p>The IP address of the instrument can be changed through the embedded web page.</p> <p>NOTE: Consult an IT professional to assist with IP settings.</p>
2	Open a web browser, such as Internet Explorer.
3	<p>In the address bar, type the current IP Address of the instrument. Press Enter.</p> <p>NOTE: The factory default IP address is set at 10.1.1.101.</p> 
4	<p>The instrument's embedded web page will display. Click the IP Configuration tab. The current IP address for the instrument is displayed. Click Make Changes.</p> 
5	<p>When prompted for a User name and Password, type ipconfig for both. Click OK.</p> 

Table 3-1 Changing the Instrument IP Address (continued)

Step	Description
6	<p>Type the desired changes to the IP address and click Submit.</p> <p>NOTE: Consult an IT professional to assist with IP settings.</p>
7	<p>A window will prompt the user to restart the instrument: Click Restart. Allow three minutes for the instrument to update configuration files.</p> <div data-bbox="277 525 349 588" data-label="Image"> </div> <div data-bbox="367 533 531 569" data-label="Section-Header"> <h3>CAUTION</h3> </div> <div data-bbox="391 602 1192 730" data-label="Text"> <p>Do not turn off the power without waiting three minutes. 3000 Micro GC must update configuration files before power is turned off. If power is turned off too soon, the files could be corrupted and render 3000 Micro GC unusable.</p> </div> <div data-bbox="623 777 1370 1285" data-label="Image"> </div>
8	<p>Connect to the instrument using the new IP address. Click (Windows) Start >> Run.</p> <div data-bbox="1169 1400 1370 1491" data-label="Image"> </div>
9	<p>In the Run window, in the Open: box, type the new IP address. Click OK.</p> <div data-bbox="826 1556 1370 1848" data-label="Image"> </div>

Table 3-1 Changing the Instrument IP Address (continued)

Step	Description
10	<p>Verify from the Command Prompt window that the PC is receiving data packets from the instrument.</p> <p>NOTE: If the ping command prompt displays Request Timed Out or Detection Limit Unreachable, then connect to the static IP address (192.168.1.99) by following the procedure in 3000 Micro GC Guide 2, Connecting to 3000 Micro GC via the Static IP Address.</p> <p>If connecting to the static IP address of the instrument is unsuccessful, contact the network administrator to check whether there is a conflict of the newly assigned IP address or any possible network issue between the instrument and the PC. Once the network issue is resolved, perform steps in this section again to configure the instrument to the new IP address.</p> <p>If connecting to the static IP address is unsuccessful, contact INFICON.</p> 

This page is intentionally blank.

3000 Micro GC Guide 4

Changing the Carrier Gas Configuration

Table 4-1 Changing the carrier gas configuration

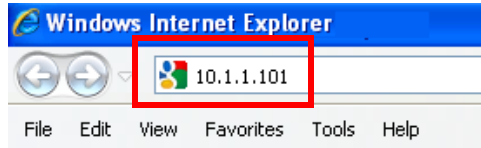
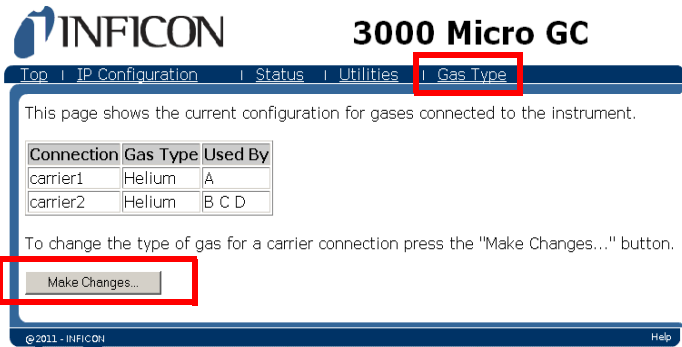

Step	Description
1	<p>The type of carrier gas used for analysis can be changed through the embedded web page.</p> <p>NOTE: If a carrier gas type is changed without following this procedure, inverted peaks will appear in the chromatogram.</p>
2	Open a web browser, such as Internet Explorer.
3	<p>In the address bar, type in the current 3000 Micro GC IP Address. Press Enter.</p> <p>NOTE: At the factory, the instrument IP address was set to 10.1.1.101.</p> 
4	<p>The instrument's embedded webpage will display. Click the Gas Type tab. The current gas configuration for installed GC modules is shown. Click Make Changes.</p> 
5	<p>When prompted for a User name and Password, type gasconfig for both.</p> 

Table 4-1 Changing the carrier gas configuration (continued)



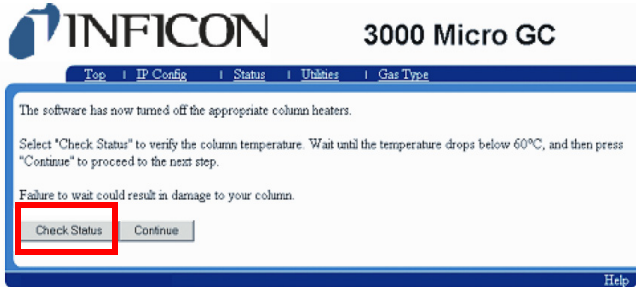
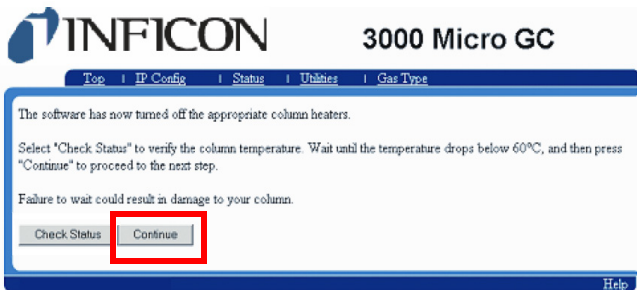

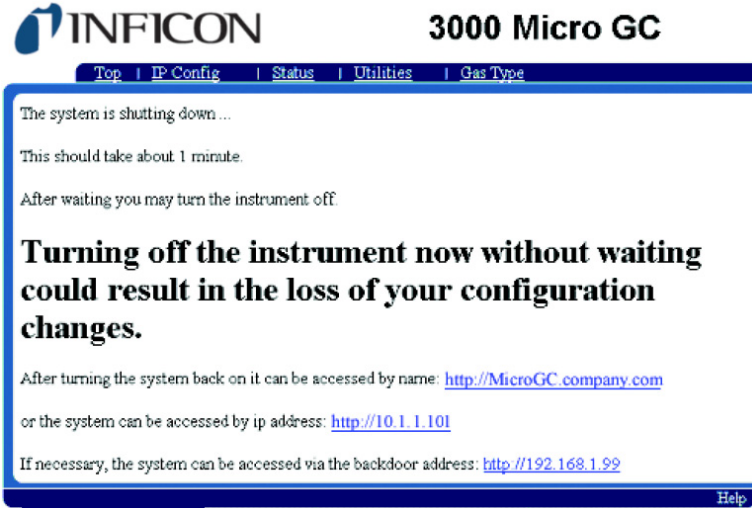
Step	Description
6	<p>Under Connection, select the carrier gas connection that corresponds to the desired GC module. Click Submit.</p> <p>NOTE: Entries m1 and m2 represent channels A and B, respectively. Entries m3 and m4 represent channels C and D, if installed.</p> <p>NOTE: In the example shown above, carrier1 is assigned to modules m1 and m2. Carrier2 is unused.</p> 
7	<p>Select the new carrier gas type from the drop-down list. Click Submit.</p> 
8	<p>The column heaters will turn off to avoid thermal shock and degradation. Click Check Status to view the current column temperatures.</p> 

Table 4-1 Changing the carrier gas configuration (continued)


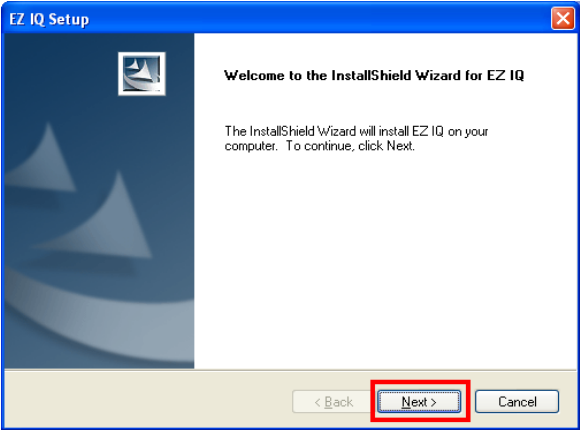
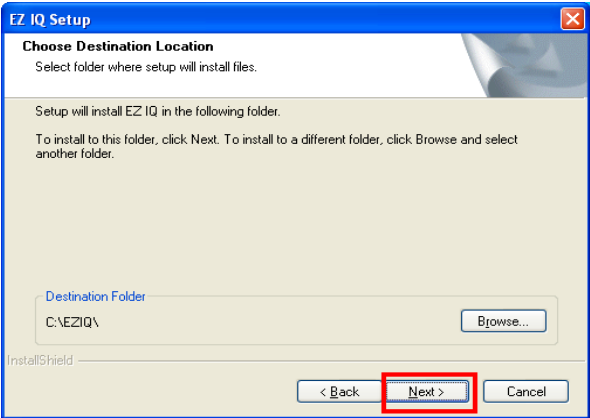
Step	Description
9	<p>When the column temperatures are below 60°C, click Continue.</p> <p>NOTE: Closing the browser or clicking back will cancel any changes.</p> 
10	<p>Change the carrier gas supply connections at the instrument back panel, and check for leaks.</p> <p>HINT: Label the new carrier gas type at the back of the instrument, for quick visual reference in the future.</p>
11	<p>A window will prompt to restart the instrument: Click Restart. Allow three minutes for the instrument to update configuration files.</p> <div data-bbox="375 961 444 1024">  </div> <p>CAUTION</p> <p>Do not turn off the power without waiting at least three minutes. 3000 Micro GC must update configuration files before power is turned off. If power is turned off too soon, the files could be corrupted and render 3000 Micro GC unusable.</p> 
12	<p>To run samples using the reconfigured gas type, click (Windows) Start >> Programs >> Chromatography >> EZ IQ Config. Update 3000 Micro GC configuration in EZ IQ. See 3000 Micro GC Guide 6, Configuring a 3000 Micro GC with EZ IQ.</p>
13	<p>A new EZ IQ method must now be created for the instrument.</p>

This page is intentionally blank.

3000 Micro GC Guide 5

Installing EZ IQ

Table 5-1 Installing EZ IQ

Step	Description
1	Insert the EZ IQ Installation CD into the PC CD-ROM drive. The CD Installer window will display. If it does not display, double-click CDInstaller.exe under the CD-ROM root drive.
2	Click Install EZ IQ to start the installation. <div></div>
3	The EZ IQ Setup - Welcome window will display. Click Next > . <div></div>
4	The EZ IQ Setup - Choose Destination Location window will display. Choose an EZ IQ software installation destination folder. The default destination folder is C:\EZIQ\ . Click Next > . <div></div>

PN 074-558-P1A

Table 5-1 Installing EZ IQ (continued)

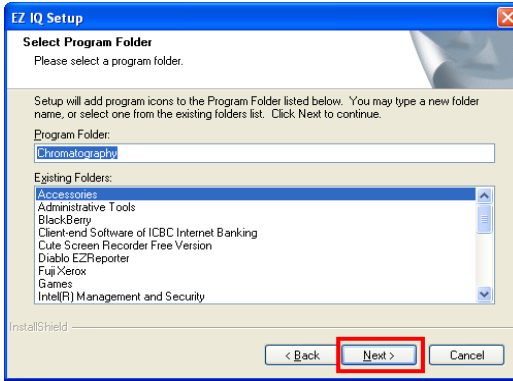
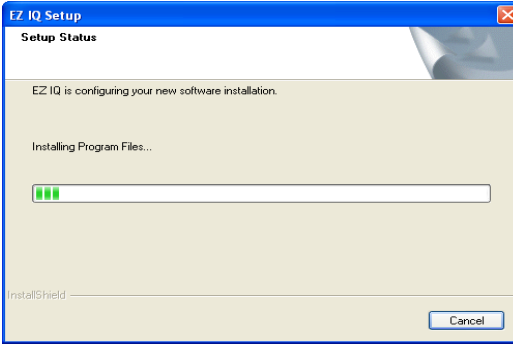
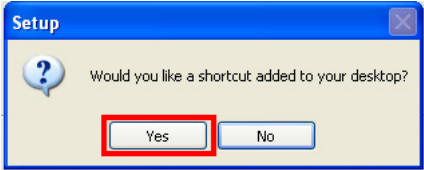
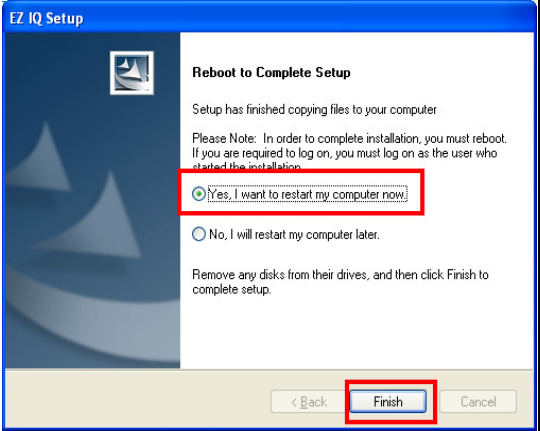
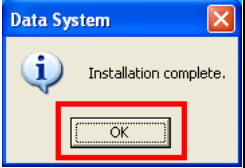
Step	Description
5	<p>The EZ IQ Setup - Select Program Folder window will display. In the Program Folder: box, type in a name for the program folder. This name will be shown in the (Windows) Start >> All Programs menu. Use the default Chromatography program folder name. Click Next >.</p> 
6	<p>The EZ IQ Setup - Setup Status window will display.</p> 
7	<p>When prompted, click Yes to create an EZ IQ program desktop short cut.</p> 

Table 5-1 Installing EZ IQ (continued)

Step	Description
8	<p>When the installation is complete, the EZ IQ Setup Reboot to Complete Setup window will display. Click Yes, I want to restart my computer now. Click Finish.</p>  The screenshot shows the 'EZ IQ Setup' window with the title 'Reboot to Complete Setup'. It contains the text: 'Setup has finished copying files to your computer. Please Note: In order to complete installation, you must reboot. If you are required to log on, you must log on as the user who started the installation.' There are two radio buttons: 'Yes, I want to restart my computer now.' (which is selected and highlighted with a red box) and 'No, I will restart my computer later.' At the bottom, there are three buttons: '< Back', 'Finish' (highlighted with a red box), and 'Cancel'.
9	<p>After the computer has restarted, the Data System window will display. Click OK to complete the installation.</p>  The screenshot shows the 'Data System' window with the title 'Data System'. It contains an information icon and the text: 'Installation complete.' Below this, there is an 'OK' button highlighted with a red box.

This page is intentionally blank.

3000 Micro GC Guide 6

Configuring a 3000 Micro GC with EZ IQ

Table 6-1 Configuring 3000 Micro GC with EZ IQ

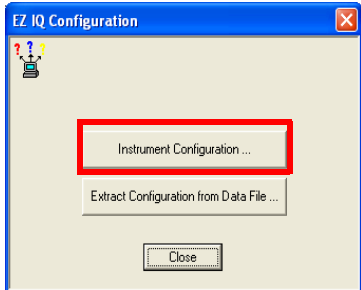
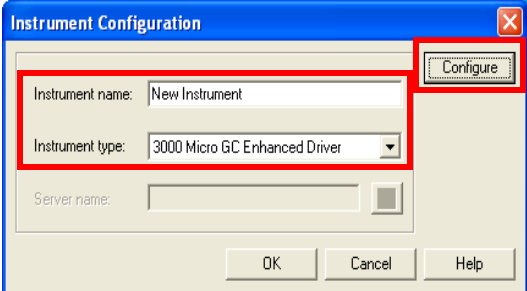
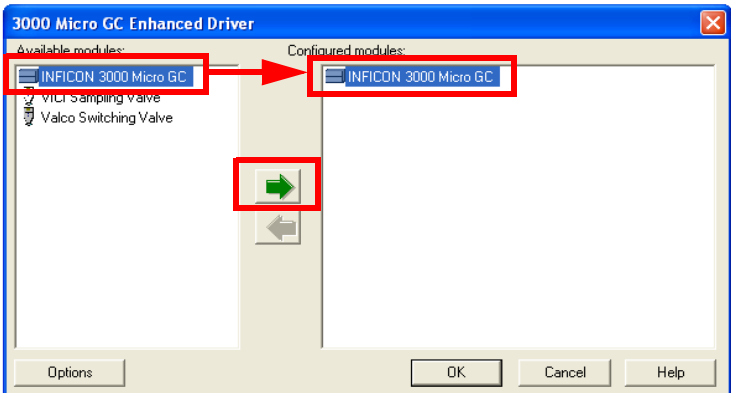
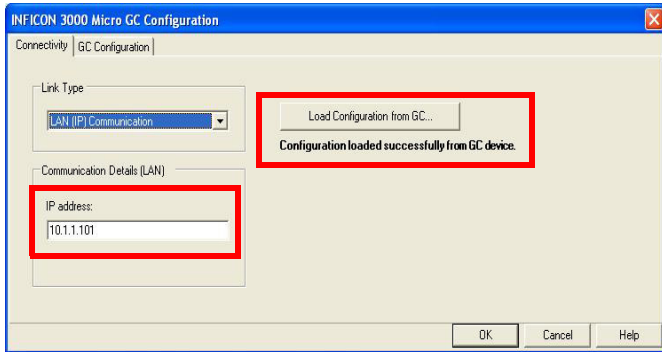
Step	Description
1	Click (Windows) Start >> All Programs >> Chromatography >> EZ IQ Config.
2	<p>The EZ IQ Configuration window will display. Click Instrument Configuration....</p> 
3	<p>On the Instrument Configuration window, type an Instrument name:, select Instrument type: 3000 Micro GC Enhanced Driver, and click Configure.</p> <p>NOTE: This guide outlines the procedure to configure the 3000 Micro GC Enhanced Driver. For the 3000 Micro GC Classic Driver, see the appropriate chapter in the PN 074-538-P1 EZ IQ Operating Manual.</p> 
4	<p>The 3000 Micro GC Enhanced Driver window will display. Under Available modules: select the INFICON 3000 Micro GC icon. Click the ➔ arrow to move the INFICON 3000 Micro GC icon to Configured modules:.</p> 

Table 6-1 Configuring 3000 Micro GC with EZ IQ (continued)

Step	Description
5	Double-click the INFICON 3000 Micro GC icon in the Configured modules panel to open the INFICON 3000 Micro GC Configuration window.
6	On the INFICON 3000 Micro GC Configuration window. Click the Connectivity tab. In the IP address: box, type the IP address of 3000 Micro GC. Click Load Configuration from GC.... Once the 3000 Micro GC configuration has been properly loaded into EZ IQ, a Configuration loaded successfully from GC device or GC Configuration Available message will display. <div data-bbox="719 535 1377 884" data-label="Image">  </div>
7	Click OK to exit the INFICON 3000 Micro GC Configuration window.
8	Click OK to exit the INFICON 3000 Micro GC Enhanced Driver window.
9	Click OK to exit the Instrument Configuration window.
10	Click Close to exit the EZ IQ Configuration window.

3000 Micro GC Guide 7

3000 Micro GC Method Setup

Table 7-1 Creating a method

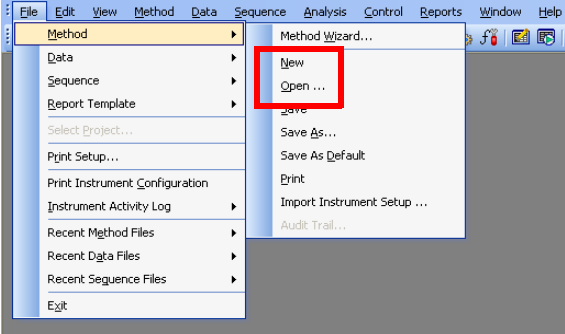
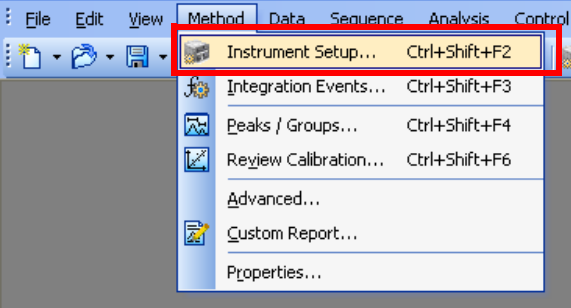
Step	Description
1	<p>It is possible to either create a new method or open an existing method in EZ IQ. Click File >> Method and select either New or Open from the drop-down list.</p> 
2	<p>To edit method parameters, click Method >> Instrument Setup.</p> 

Table 7-1 Creating a method (continued)

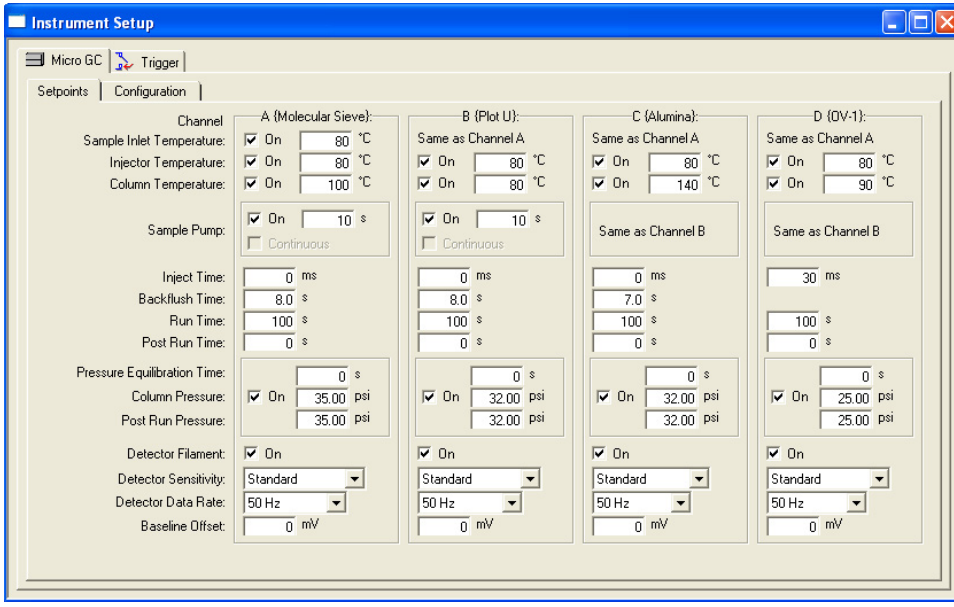
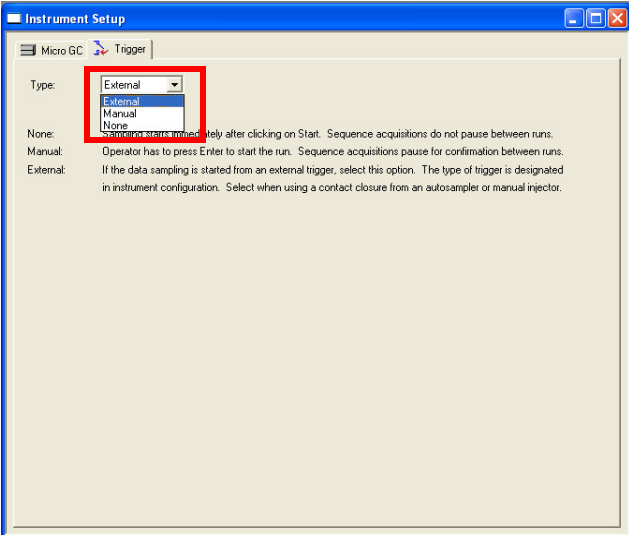
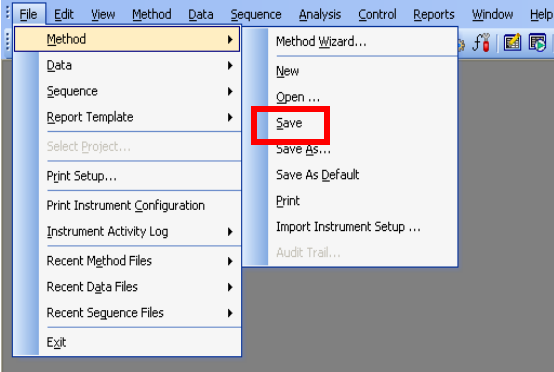
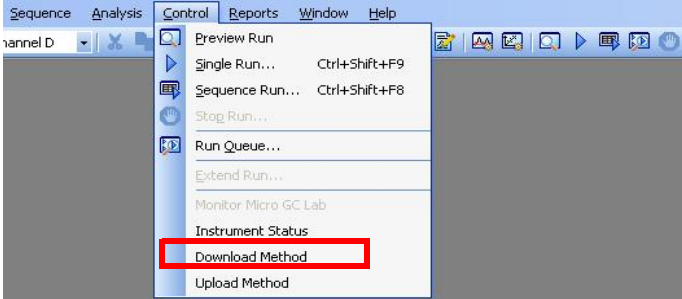
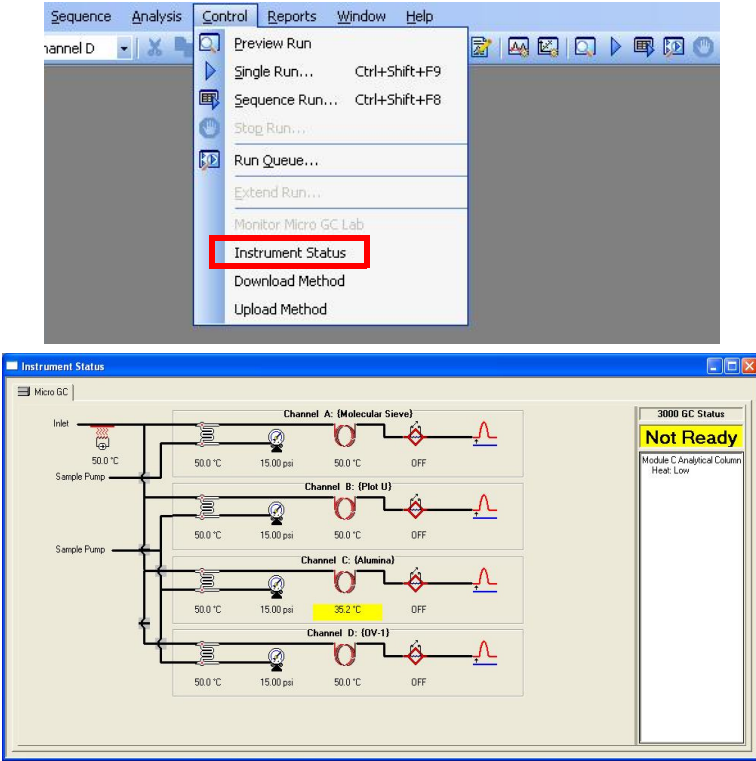
Step	Description
3	<p>The Instrument Setup window will display. It will display parameters that can be edited.</p> <p>NOTE: The Instrument Setup window is configuration dependent.</p> <p>Populate sample acquisition parameters as desired, such as Injector Temperature, Column Temperature, Sample Pump time, Column Pressure, etc. Ensure that the Run Time is long enough for the last expected peak to elute. If the elution time is unknown, conduct a run and verify the presence of the last expected peak. If necessary, increase the Run Time.</p> <p>The default Inject Time should be set between:</p> <ul style="list-style-type: none"> • 0-50 ms for a fixed volume injector • 0-100 ms for a variable volume injector • 0-350 ms for a variable large volume injector (LVI) • 0-250 ms for a backflush injector 

Table 7-1 Creating a method (continued)

Step	Description
4	<p>Click the Trigger tab. In the Type: list, select the trigger type.</p> <ul style="list-style-type: none"> If the sample run will be initiated directly through EZ IQ, select None. If the sample run will be triggered by an external event through a remote control connection, select External. 
5	<p>To save the method, click File >> Method >> Save to overwrite the existing method file or Save As to save a new method file.</p> 
6	<p>When a new method is created, or when a method is changed or opened, it must be downloaded to 3000 Micro GC. Click Control >> Download Method.</p> 

PN 074-558-P1A

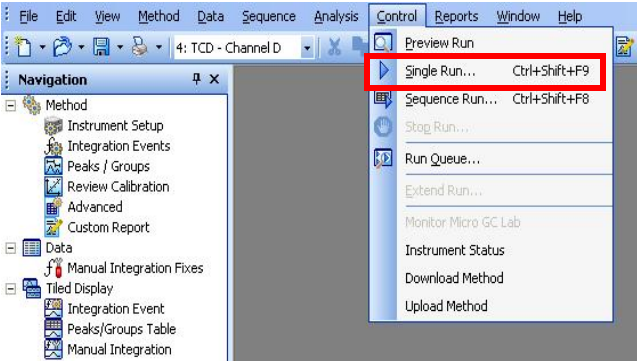
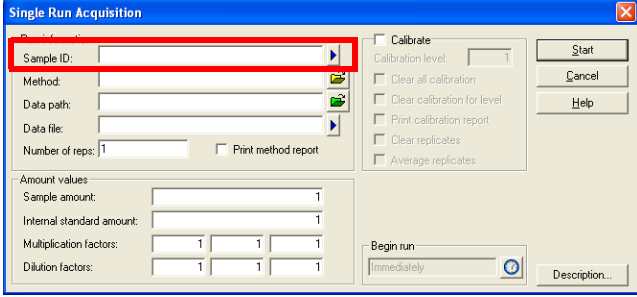
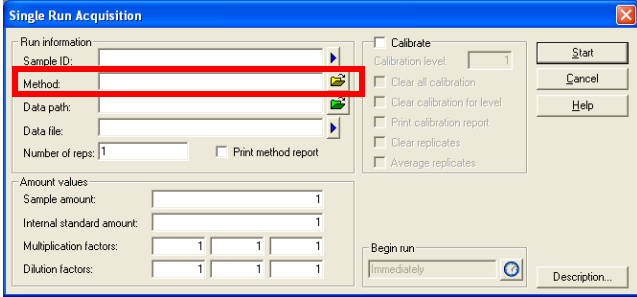
Table 7-1 Creating a method (continued)

Step	Description
7	<p>Click Control >> Instrument Status to view the status of the instrument.</p> <p>NOTE: The status will read Not Ready, highlighted in yellow, if parameters have not reached their set points. Parameters that are changing are highlighted in yellow. A run cannot be started unless the status reads Ready, highlighted in green.</p>  <p>The screenshot shows the 'Instrument Status' window with four channels listed: Channel A (Molecular Sieve), Channel B (Plot U), Channel C (Alumina), and Channel D (OV-1). Channel C is highlighted in yellow, indicating it is not ready. The status bar on the right shows 'Not Ready' in yellow. The parameters for Channel C are: 50.0 °C, 15.00 psi, 35.2 °C (highlighted in yellow), and OFF. The status bar also shows 'Module C Analytical Column Heat Low'.</p>

3000 Micro GC Guide 8

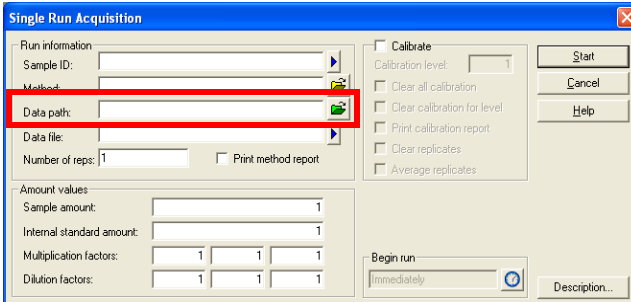
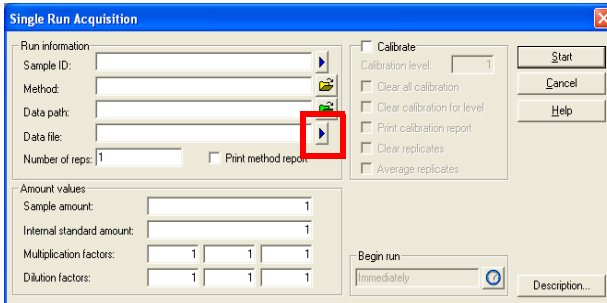
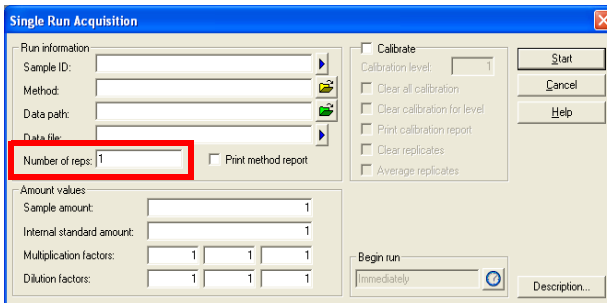
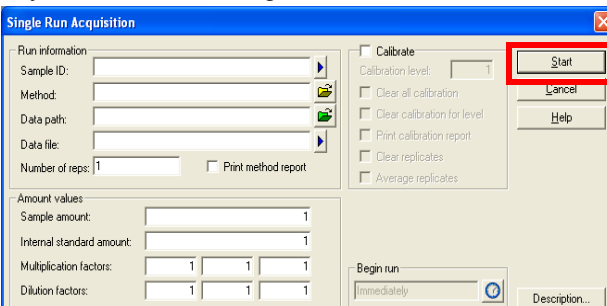
Running a Sample

Table 8-1 Running a sample

Step	Description
1	<p>To start the run, click Control >> Single Run.</p> 
2	<p>The Single Run Acquisition window will display. In the Sample ID: box, type a Sample ID.</p> 
3	<p>The method will default to the method that is currently loaded. To change the method, click the folder icon located on the right of the Method: box and select the desired method.</p> 

PN 074-558-P1A

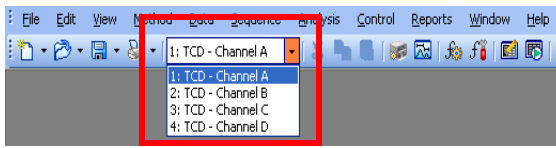
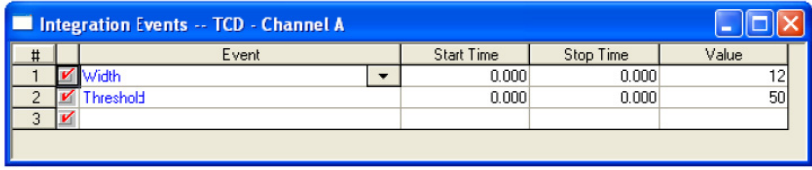
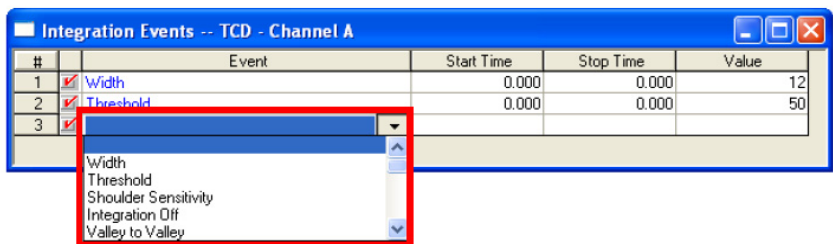
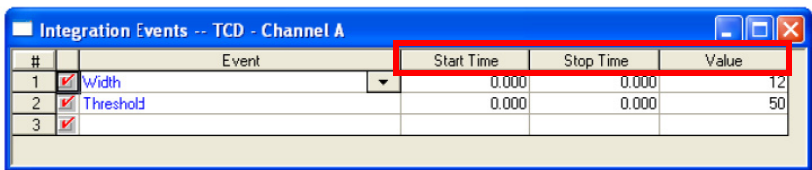
Table 8-1 Running a sample (continued)

Step	Description
4	<p>Select a path for storage of data files by selecting the folder icon located on the right of the Data path: box.</p> 
5	<p>The Data file: box is used to create the filename. To automatically include certain descriptors such as Sample ID, Method Name, and Date and Time, click the blue arrow, located to the right of the Data File field. Select the desired descriptors.</p> 
6	<p>In the Number of Reps box, type the desired number of repetitions.</p> 
7	<p>Click Start to begin a run. Data will be displayed in the chromatogram window.</p> 

3000 Micro GC Guide 9

Integrating Chromatograms

Table 9-1 Integrating Chromatograms


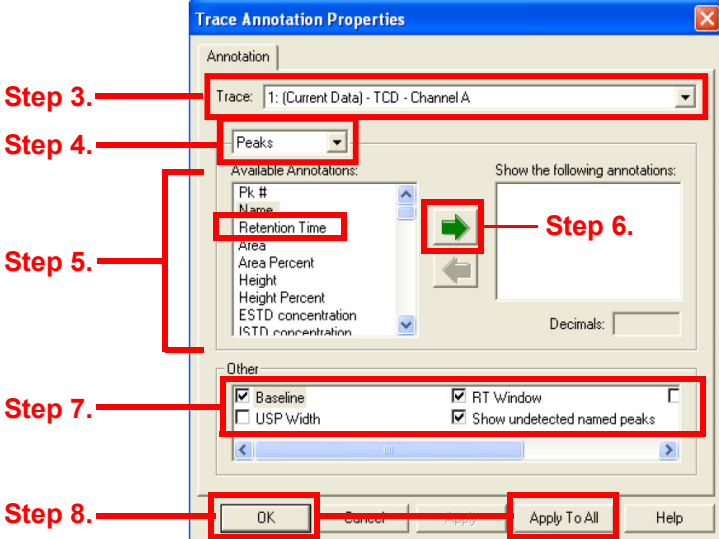
Step	Description
1	Integration events are used to customize the integration of certain peaks or regions of the chromatogram. Refer to the 074-538 EZ IQ Operating Manual or online help in the EZ IQ software for descriptions of individual events. In EZ IQ , click Method >> Integration Events .
2	Select a channel to add or edit integration events. To change the channel, select the appropriate channel from the drop-down list on the toolbar. 
3	The Integration Events window will display the default integration events. 
4	To add a new event, click on a blank field under Event in the table. Select the desired integration event from the drop-down list. 
5	Adjust the integration Start Time , Stop Time , and Value as desired. 
6	Click Analysis >> Analyze to apply the changes.
7	To save the method, click File >> Method >> Save .

This page is intentionally blank.

3000 Micro GC Guide 10

Setting Annotations

Table 10-1 Setting Annotations

Step	Description
1	<div>In EZ IQ, right-click on a chromatogram and click Annotations.</div> <div></div>
2	<div>The Trace Annotation Properties window will appear.</div> <div></div>
3	In the Trace: list, select the desired chromatogram.
4	For the selected chromatogram, select Peaks or Groups from the drop-down list.
5	On the Available Annotations: list, select the features to annotate. For example, select Retention Time to display the retention time above the integrated peaks.
6	<div>When an annotation is highlighted, double-click the annotation or click the green arrow key to move it from Available Annotations: to Show the following annotations:.</div> <div>NOTE: For certain annotations, the number of decimal places can be designated. Enter this value in the Decimals box for the highlighted item.</div>

PN 074-558-P1A

Table 10-1 Setting Annotations (continued)

Step	Description
7	Select the check boxes to display Baseline , USP Width , Retention Time Window , Show Undetected Named Peaks , or Group Range on the chromatogram.
8	When finished selecting annotations, click Apply to All or OK .

3000 Micro GC Guide 11

Qualifying Peaks and Groups

Table 11-1 Qualify peaks and groups

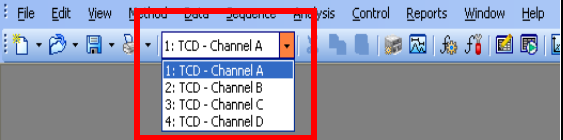
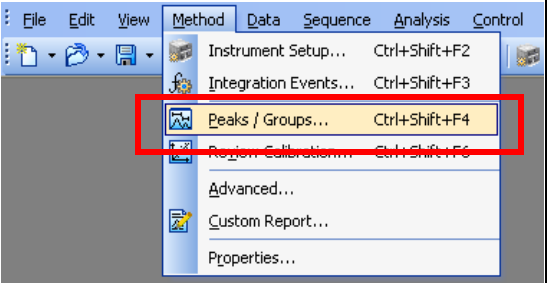
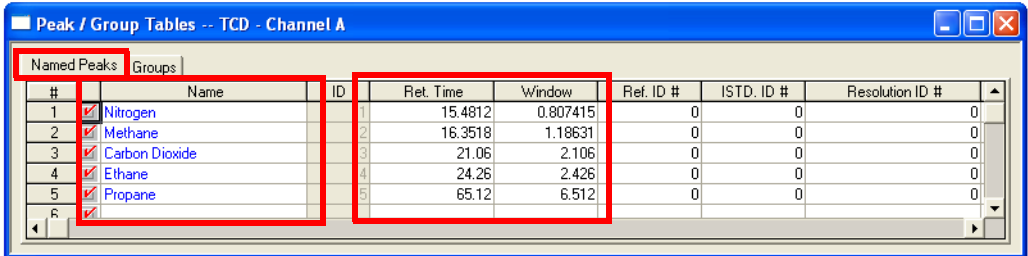
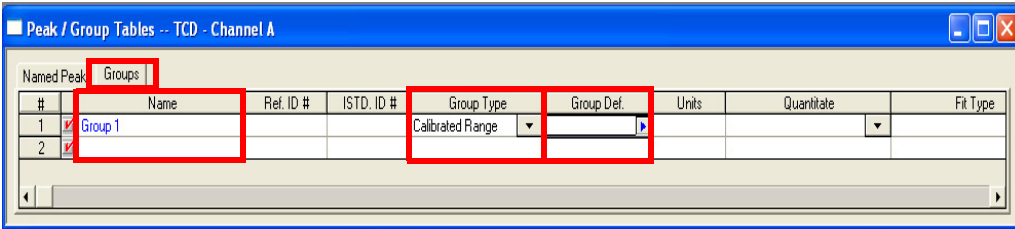
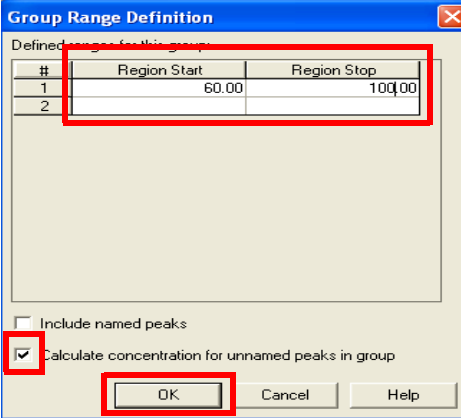
Step	Description
1	<p>Once a run is finished, peaks and groups can be identified based on retention time. Select a channel to add or edit peak or group information from the drop-down list on the tool bar.</p> 
2	<p>Click Method >> Peaks/Groups.</p> 

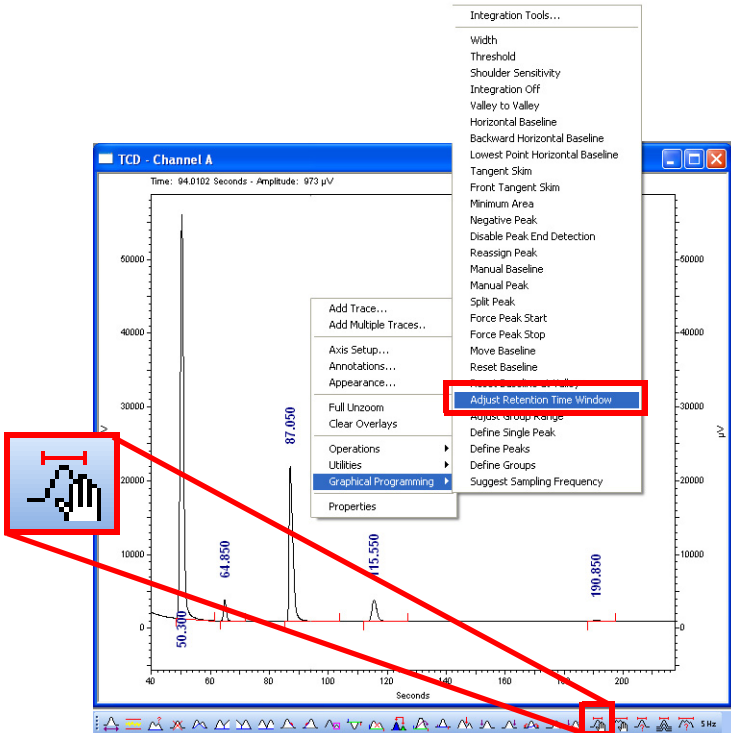
Table 11-1 Qualify peaks and groups (continued)

Step	Description
3	<p>The Peak/Group Tables window will display. There are tabs for Named Peaks and Groups. On the Named Peaks tab, enter the Name, Ret.Time, and retention time Window for each component in the channel.</p>  <p>To enter group information, click the Groups tab. In a blank Name box, type the name of the group. In the Group Type list, select Calibrated Range.</p>  <p>Click the blue arrow on the right of the Group Def. field. The Group Range Definition window will display. Type a Region Start and Region Stop, in minutes or seconds.</p> <p>NOTE: In the example, Region Start and Region Stop are displayed in seconds.</p> <p>Ensure that Calculate Concentrations for Unnamed Peaks in Group is selected from the drop-down list. Click OK to return to the Groups tab.</p> 
4	Repeat Steps 1 to 3 for all channels.
5	Click Analysis >> Analyze to apply the changes.
6	Save the method by clicking File >> Method >> Save .

3000 Micro GC Guide 12

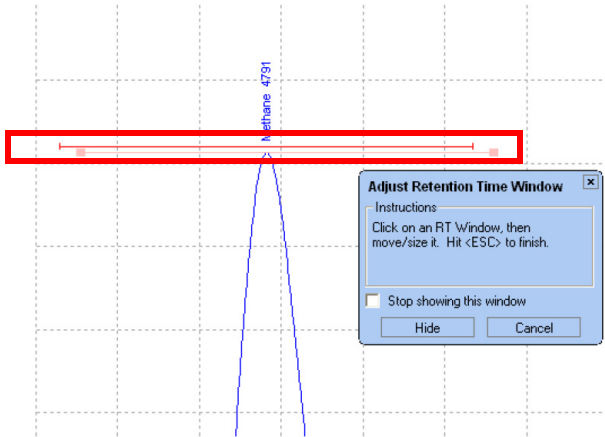
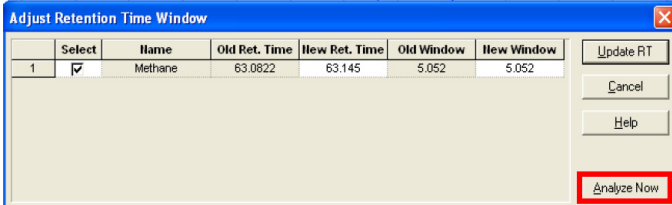
Adjusting Retention Time Windows

Table 12-1 Adjusting retention time windows

Step	Description
1	<p>In EZ IQ, click the Adjust Retention Time Window icon on the Int. Events tool bar. Alternatively, right-click on the chromatogram and click Graphical Programming >> Adjust Retention Time Window.</p> 

PN 074-558-P1A

Table 12-1 Adjusting retention time windows (continued)

Step	Description
2	<p>A blue Adjust Retention Time Window window will display. Click directly on the desired retention time window.</p> <p>HINT: Click and drag from the corner of the window to resize.</p> 
3	<p>Press <Esc> to display the Adjust Retention Time Window option. Click Analyze Now to apply the changes and exit the window.</p> 

3000 Micro GC Guide 13

Conducting a Single Level Calibration

Table 13-1 Conduct a single level calibration

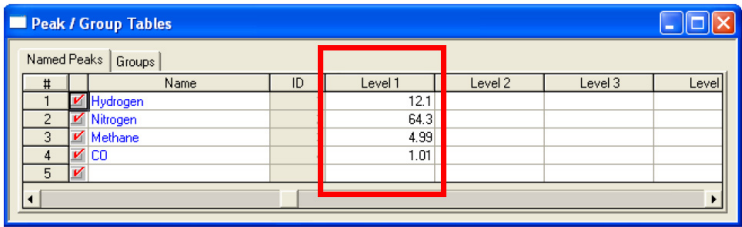
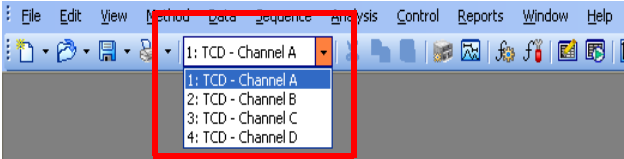
Step	Description
1	<p>Prior to conducting qualitative or quantitative analysis on 3000 Micro GC, it is necessary to properly calibrate the instrument using a known calibration gas. The calibration process identifies the elution order of the component peaks in a chromatogram, as well as the detector response each component demonstrates with the instrument.</p> <ul style="list-style-type: none"> If the sample gas concentration is close to the calibration gas, a single level calibration is sufficient. Go to step 2. If the sample gas concentration varies over a broad range, a multiple level calibration is recommended. See 3000 Micro GC Guide 14, Conducting a Multiple Level Calibration.
2	To conduct a single level calibration, run a sample of a calibration gas (i.e., a sample with known component concentrations). The generated data file can be used to define the calibration peaks.
3	Connect the calibration gas to the front inlet of 3000 Micro GC. Ensure that the calibration gas output pressure and the 3000 Micro GC inlet is properly regulated to between 1-25 PSI. Open the gas cylinder to begin flow to the instrument.
4	Click Control >> Single Run to run the sample with the currently loaded method. Type the desired Sample ID and click Start . After the run is complete, ensure all peaks are integrated and annotated.
5	<p>The known concentration of components in the calibration gas must be recorded in the Peaks/Groups table.</p> <p>Click Method >> Peaks/Groups. The Peak/Group Tables window will display. Use the scroll bar to scroll until the Level 1 column is visible.</p> 
6	<p>Enter the concentration for the component under Level 1. Repeat this for all peaks and groups in each of the channels.</p> <p>To add or edit calibration information, select a channel from the drop-down list on the tool bar.</p> 

Table 13-1 Conduct a single level calibration (continued)

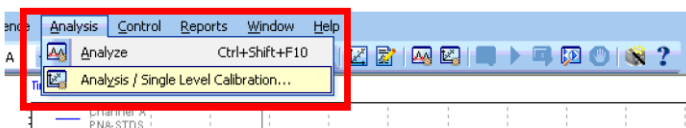
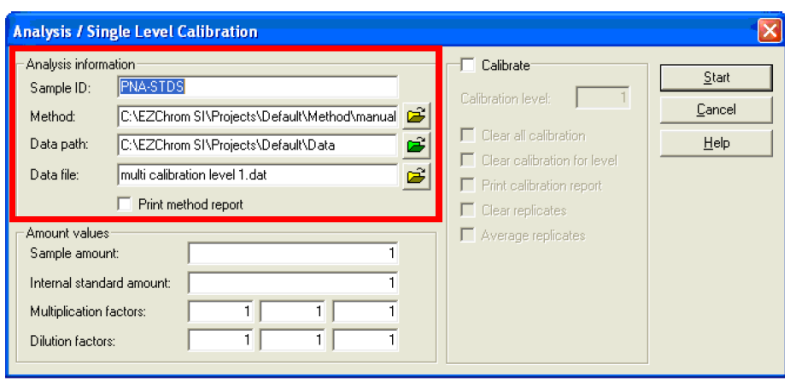
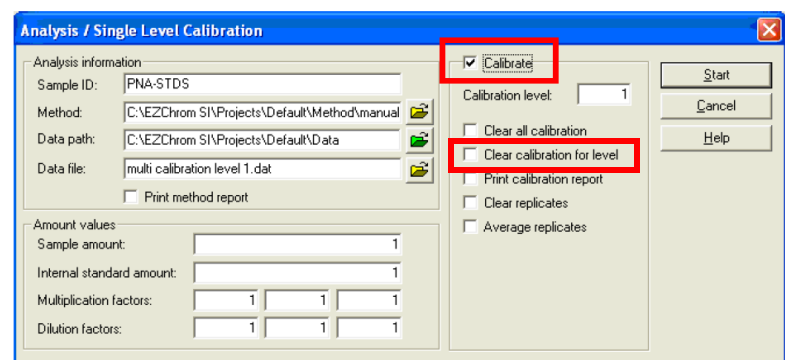
Step	Description
7	To calibrate, click Analysis >> Analysis/Single Level Calibration . 
8	The Analysis/Single Level Calibration window will display. 
9	In the Sample ID: box, type the sample identification, if not already present.
10	In the Method: box, select the name of the method to be calibrated, if not already present. To do so, click the Open file icon adjacent to the field.
11	In the Data path: box, select the data path name, if not already present. To do so, click the Open file icon adjacent to the field.
12	In the Data file: box, select the file name, if not already present. To do so, click the Open file icon adjacent to the field.
13	The Amount values will remain set at 1 ; a value of 1 corresponds to a single level calibration.
14	Select Calibrate . In the Calibration Level box, type 1 . NOTE: For the initial calibration, it is not necessary to select any of the boxes involving clearing calibrations or replicates. However, if the method was previously calibrated and the calibrated information needs to be replaced, select Clear all calibration before starting. 

Table 13-1 Conduct a single level calibration (continued)

Step	Description
15	When finished, click Start .
16	When the analysis is complete, the chromatogram will be integrated, and the areas for the peaks identified as calibration compounds will be entered into the method. In addition, a calibration curve will be generated using these areas. The method is now calibrated for a single level.
17	To save the calibration to the method, click File >> Method >> Save .

This page is intentionally blank.

3000 Micro GC Guide 14

Conducting a Multiple Level Calibration

Table 14-1 Conduct a multiple level calibration

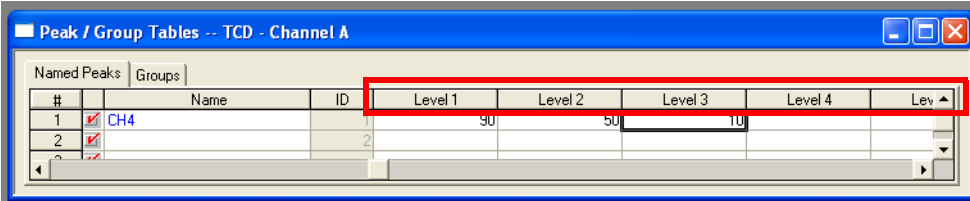
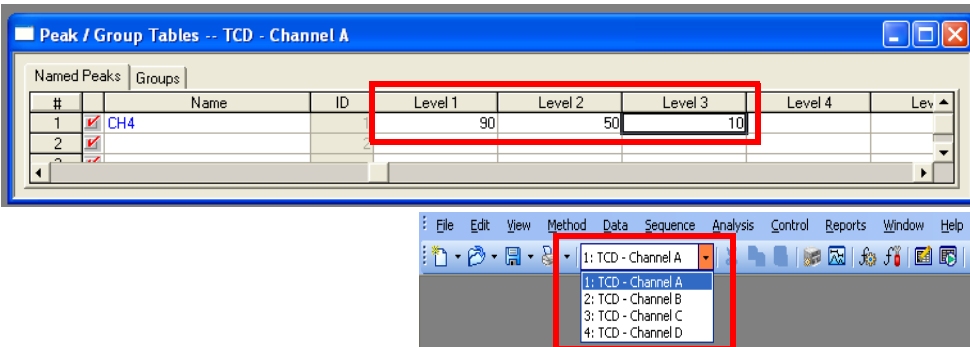
Step	Description
1	If a multiple level calibration is required, run sample calibration standards of different concentration levels.
2	<p>The known concentration of components in the calibration gas must be recorded in the Peaks/Groups table.</p> <p>Click Method >> Peaks/Groups. The Peak/Group Tables window will display. Use the scroll bar to display the columns for Level 1, Level 2, etc.</p> 
3	<p>Type the concentrations for the component at each level. Repeat this for all peaks and groups in each of the channels.</p> <p>NOTE: Each level represents one calibration gas standard.</p> <p>Select a channel to add or edit calibration information via the dropdown list on the tool bar.</p> 
4	<p>Connect the calibration gas. Ensure that the tank is properly regulated to between 1-25 PSI, if going direct to the 3000 Micro GC inlet. Turn on the desired calibration gas.</p> <p>Connect the calibration gas to the front inlet of 3000 Micro GC. Ensure that the calibration gas output pressure and the 3000 Micro GC inlet is properly regulated to between 1-25 PSI. Open the gas cylinder to begin gas flow to the instrument.</p>
5	Click Control >> Single Run to run the sample on the desired method. Type the Sample ID into the box and click Start . After the run is complete, ensure all peaks are integrated and annotated.

Table 14-1 Conduct a multiple level calibration (continued)

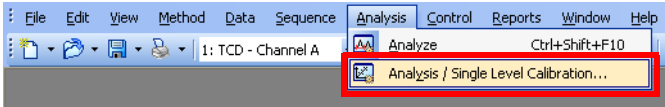
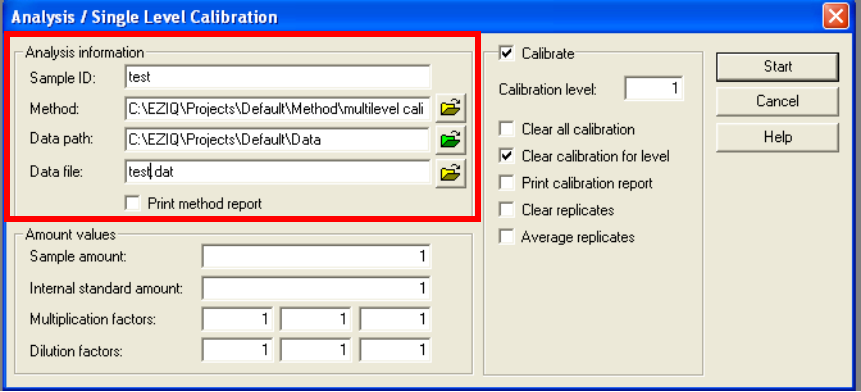
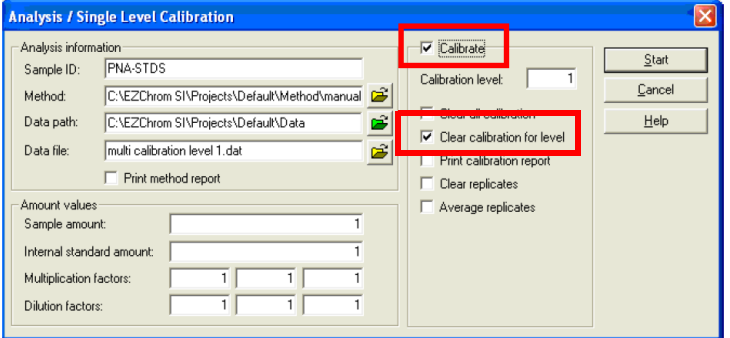
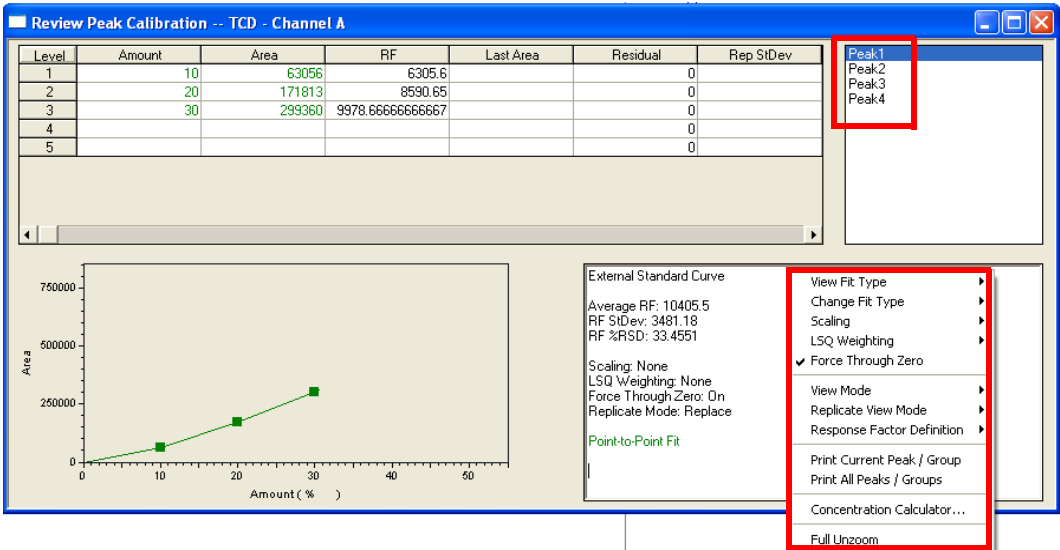
Step	Description
6	To calibrate, click Analysis >> Analysis/Single Level Calibration . 
7	The Analysis/Single Level Calibration window will display. 
8	In the Sample ID : box, type the sample identification, if not already present.
9	In the Method : box, select the name of the method to be calibrated, if not already present, by clicking the Open file icon adjacent to the field.
10	In the Data path : box, select the data path name, if not already present, by clicking the Open file icon adjacent to the field and selecting the desired data path.
11	In the Data file : box, select the file name, if not already present, by clicking the Open file icon adjacent to the field.
12	The Amount values are to remain set to 1 ; a value of 1 corresponds to a single level calibration.
13	Select Calibrate . In the Calibration level : box, type 1 . NOTE: For the initial calibration, it is not necessary to select any of the boxes involving clearing calibrations or replicates. However, if the method was previously calibrated and the calibrated information needs to be replaced, select Clear calibration for level before starting. 
14	Click Start after completing the window.

Table 14-1 Conduct a multiple level calibration (continued)

Step	Description
15	When the analysis is complete, the chromatogram will be integrated. The areas for the peaks identified as calibration compounds will be entered into the method. In addition, a calibration curve will be generated using these areas. The method is now calibrated for one of the desired levels.
16	Repeat steps 4 to 15 to calibrate other levels, one level at a time.
17	After completing a calibration in the window, click Method >> Review Calibration to review the calibration curve and its associated data. Each peak will have its own calibration curve. If desired, right-click on a blank area to bring up options such as View Fit Type , Change Fit Type , and a Concentration Calculator .



This page is intentionally blank.

3000 Micro GC Guide 15

Calibration with Average Replicates

Table 15-1 Conducting a calibration with replicates

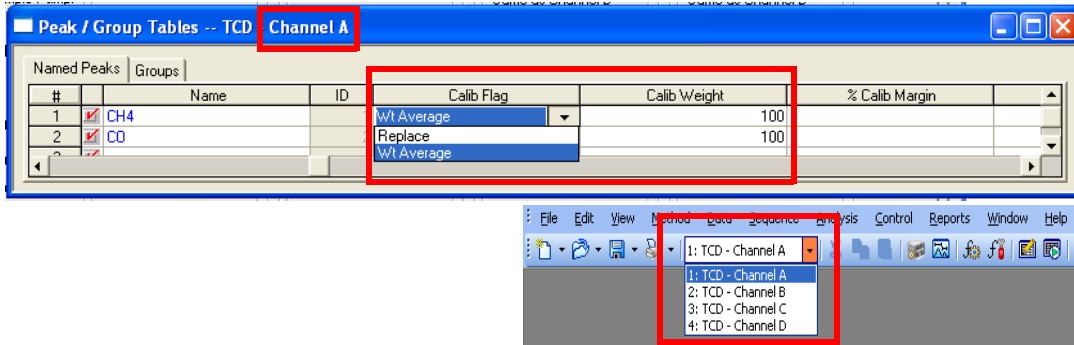
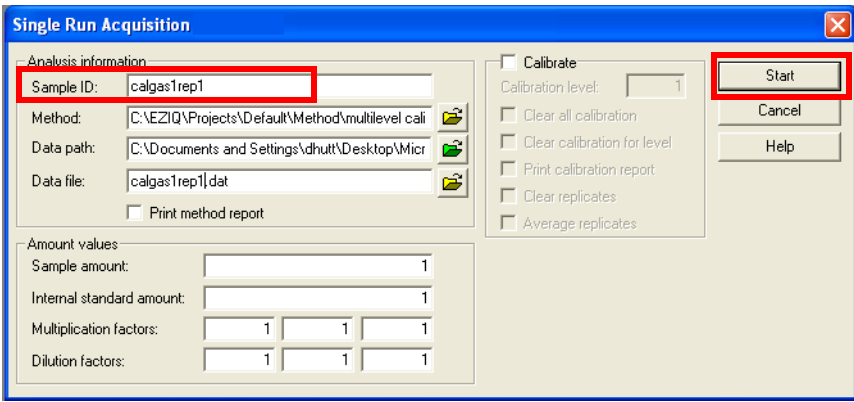
Step	Description
1	This guide will go through the procedure for a three point calibration of a single level. The procedure can be repeated for multiple levels, if desired.
2	<p>Click Method >> Peaks/Groups. The Peak/Group Tables window will display. Use the scroll bar to display the Calib Flag and Calib Weight columns.</p> <p>Change the Calib Flag to Wt Average using the drop-down list.</p> <p>Enter 100 for the Calib Weight. Repeat for each component in each channel.</p> <p>NOTE: Select a channel from the drop-down list on the toolbar to add or edit calibration information.</p> 
3	Connect the calibration gas to the front inlet of 3000 Micro GC. Ensure that the calibration gas output pressure is properly regulated between 1-25 PSI. Open the gas cylinder to begin gas flow to the instrument.
4	<p>Click Control >> Single Run. In the Sample ID: box, type a unique sample identifier. In this example, the Sample ID: is calgas1rep1. Click Start to begin a run.</p> 
5	After the run is complete, ensure all peaks are integrated and annotated.

Table 15-1 Conducting a calibration with replicates (continued)

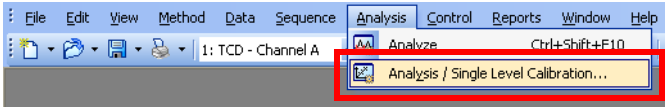
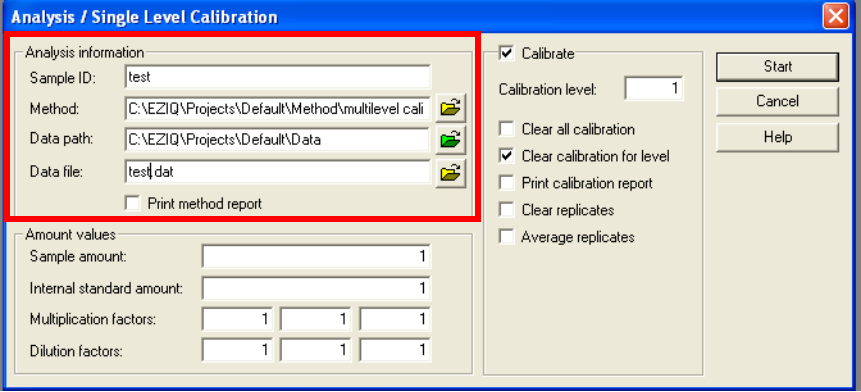
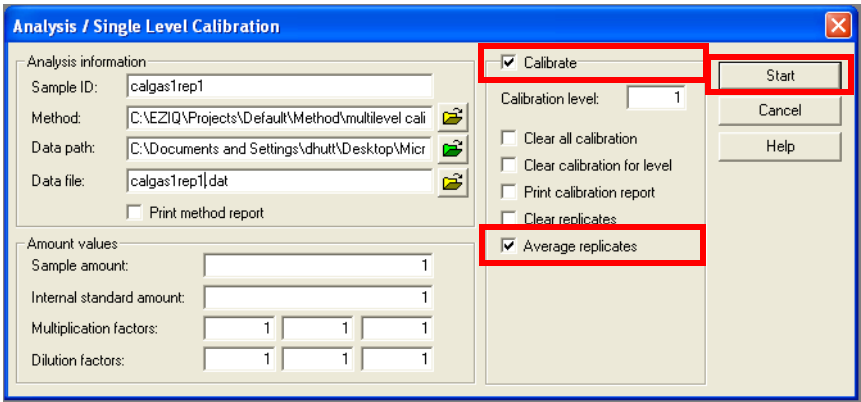
Step	Description
6	To calibrate, click Analysis >> Analysis/Single Level Calibration . 
7	The Analysis/Single Level Calibration window will display. 
8	In the Sample ID: box, type the sample identification, if not already present.
9	In the Method: box, select the name of the method to be calibrated, if not already present, by clicking the Open file icon adjacent to the field.
10	In the Data path: box, select the data path name, if not already present, by clicking the Open file icon adjacent to the field and selecting the desired data path.
11	In the Data file: box, select the file name, if not already present, by clicking the Open file icon adjacent to the field.
12	The Amount values are to remain set to 1 ; a value of 1 corresponds to a single level calibration.
13	Select the Calibrate setting and type 1 for Calibration Level . Select Average replicates . NOTE: For the initial calibration, it is not necessary to select boxes that involve clearing calibrations or replicates. However, if the method was previously calibrated and the calibrated information needs to be replaced, select Clear all calibration . 

Table 15-1 Conducting a calibration with replicates (continued)

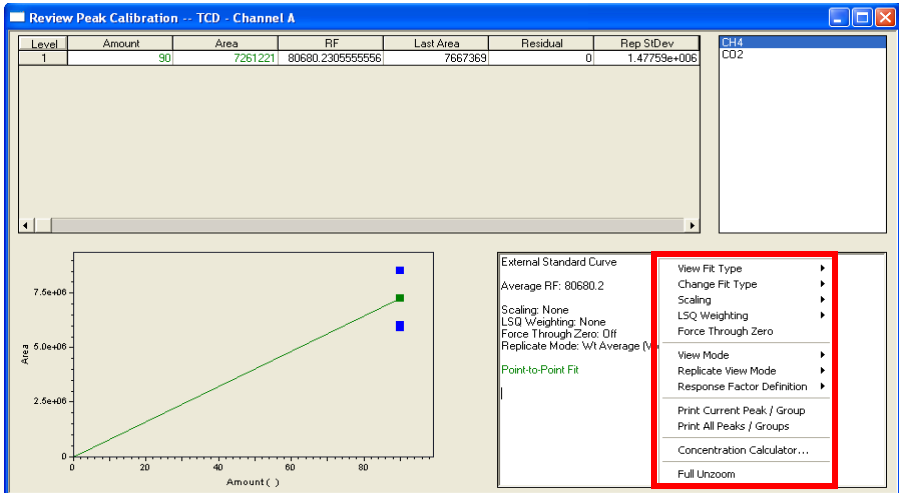
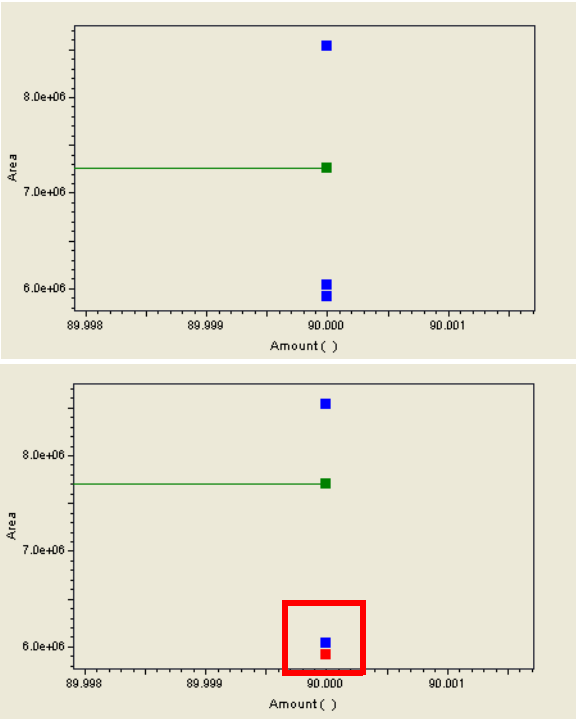
Step	Description
14	When finished, click Start .
15	When the analysis is complete, the chromatogram will be integrated, and the areas for the peaks identified as calibration compounds will be entered into the method.
16	With the calibration gas still connected and flowing, repeat steps 4 through 14 for each replicate. Ensure that each Sample ID is unique, for example calgas1rep2 , calgas1rep3 , etc.
17	<p>After completing a calibration in the window, a calibration curve and associated data can be reviewed by clicking Method >> Review Calibration. Each peak will have its own calibration curve. If desired, right click on a black area on the window and select commands such as View Fit Type, Change Fit Type, and a Concentration Calculator.</p> <p>The peak Area will be reported as the average area of the replicates.</p> 

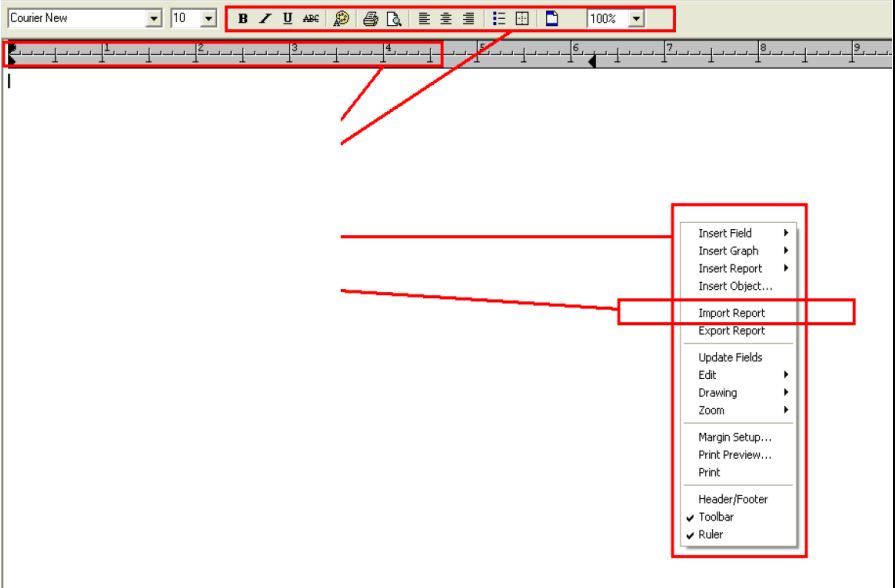
Table 15-1 Conducting a calibration with replicates (continued)

Step	Description
18	<p>On the calibration curve, a blue point represents a single replicate, and a green point represents the average of the replicates. Left-clicking on a single blue point will turn the point red and remove it from the average calculation.</p> 

3000 Micro GC Guide 16

Customizing a Method Report

Table 16-1 Customize a report

Step	Description
1	The default method report is blank. To add fields, graphs, and data, click Method >> Custom Report .
2	A method report window will display. 
3	Sections can be added to the report by right-clicking anywhere on the method report. Select the desired options such as Insert Field , Insert Graph , Insert Report , or Insert Object .
4	EZ IQ contains various pre-made template files. To import a template file, right-click in the method report window. Select Import Report .

PN 074-558-P1A

Table 16-1 Customize a report (continued)

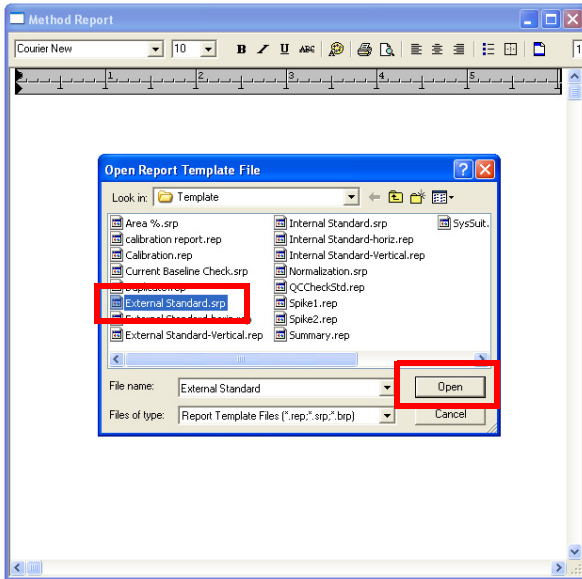
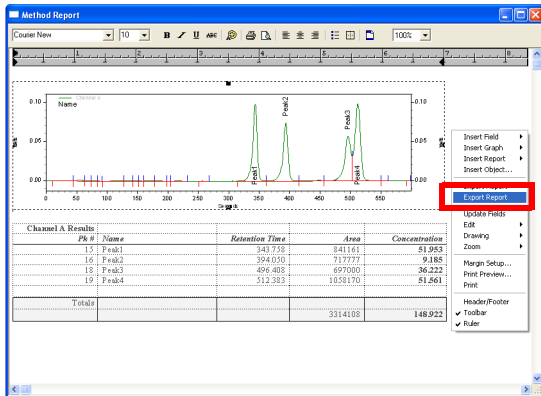
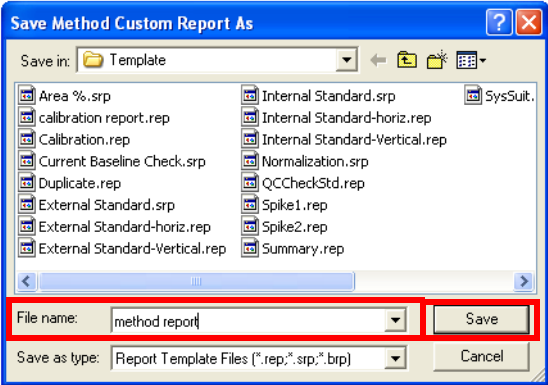
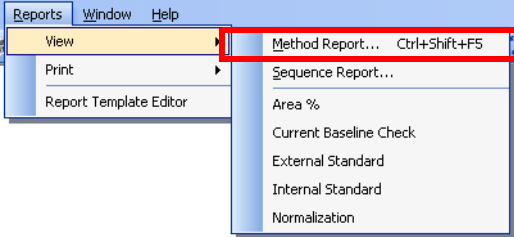
Step	Description
5	<p>The Open Report Template File window will prompt to open a template file, such as External Standard.srp, shown below. Click Open to load the template file.</p> 
6	<p>To edit the template, right-click and insert options such as Insert Field, Insert Graph, Insert Report, and Insert Object.</p>
7	<p>Once all desired information has been modified or added, right-click in the Method Report and select Export Report.</p> 

Table 16-1 Customize a report (continued)

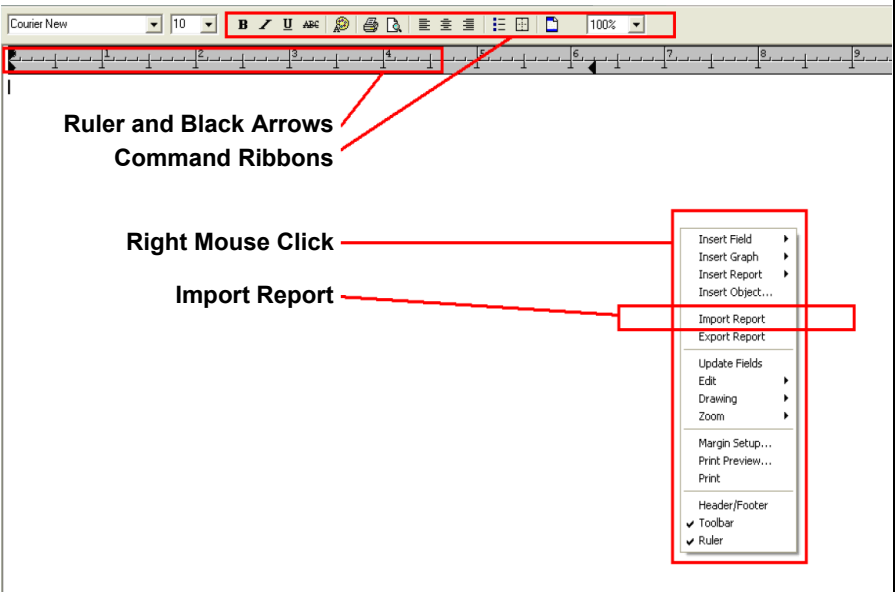
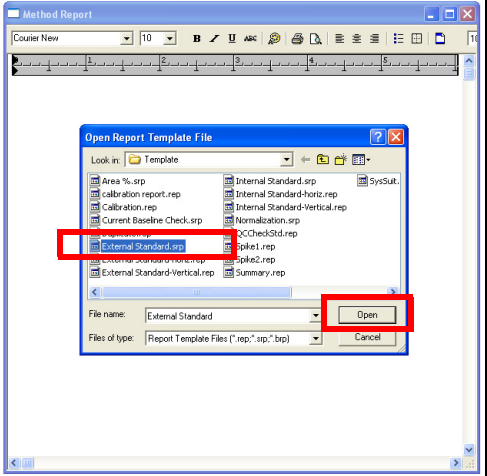
Step	Description
8	<p>The Save Method Custom Report As window will display. In the File Name: box, type a method report name. Click Save.</p> 
9	<p>To view the custom method report, select Reports >> View >> Method Report.</p> 
10	<p>To print the custom Method Report, right-click in the Method Report window, then select Print from the drop-down list.</p>

This page is intentionally blank.

3000 Micro GC Guide 17

Customizing an ESTD Report

Table 17-1 Customizing an external standard (ESTD) report

Step	Description
1	An External Standard (ESTD) Report contains information on retention time, area, and mole percent concentration. It is not method-specific. The same template will be applied to any method that is opened.
2	Click Method >> Custom Report . A method report window will display. <div></div>
3	To import the ESTD Report file, right-click in the method report window, and select Import Report from the dropdown list.
4	The Open Report Template File window will display. Select External Standard.srp and click Open to load the template file. <div></div>

PN 074-558-P1A

Table 17-1 Customizing an external standard (ESTD) report (continued)

Step	Description
5	To edit the report, right-click and select options such as Insert Field , Insert Graph , Insert Report , and Insert Object .
6	Once all desired information has been modified or added, right click in the Method Report window and select Export Report .
7	The Save Method Custom Report As window will open. Click External Standard.srp . Click Save . Click Yes to overwrite the current file.

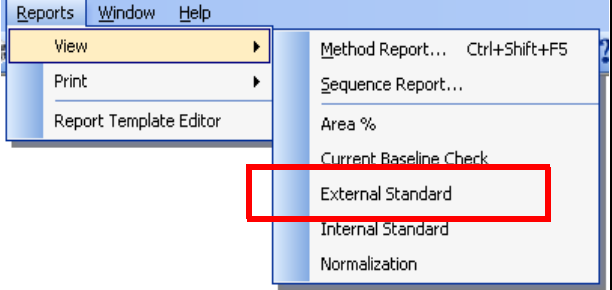
The screenshot shows the 'Method Report' window. At the top is a toolbar with various icons. Below it is a chromatogram plot with peaks labeled Peak1, Peak2, Peak3, and Peak4. Below the plot is a table titled 'Channel A Results'.

PK #	Name	Retention Time	Area	Concentration
15	Peak1	343.758	841161	51.953
16	Peak2	394.050	717777	9.185
18	Peak3	496.408	697000	36.222
19	Peak4	512.383	1058170	51.561
Totals			3314108	148.922

A right-click context menu is open on the right side of the window, with the 'Export Report' option highlighted in blue.

The screenshot shows the 'Save Method Custom Report As' dialog box. The 'Save in' field is set to 'Template'. The file name is 'External Standard'. The 'Save as type' is 'Report Template Files (*.rep;*.srp;*.trp)'. A warning message is displayed: 'C:\E2D\Projects\Default\Template\External Standard.srp already exists. Do you want to replace it?' with 'Yes' and 'No' buttons. The 'Yes' button is highlighted with a red box.

Table 17-1 Customizing an external standard (ESTD) report (continued)

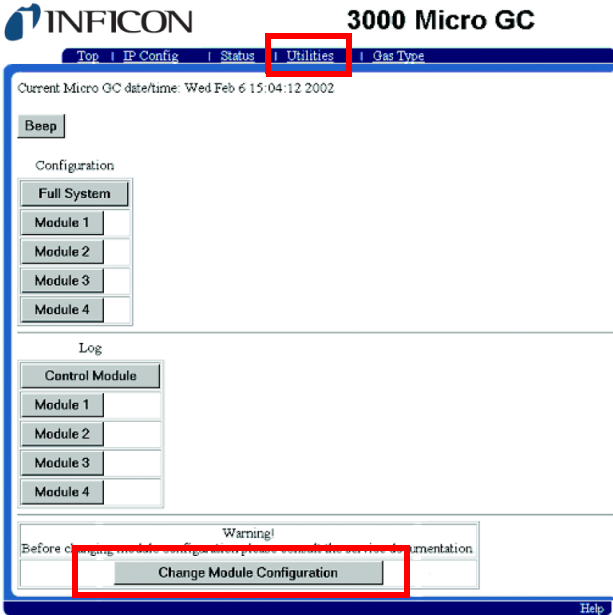
Step	Description
8	<div><div>To view the ESTD report, click Reports >> View >> External Standard.</div><div></div></div>
9	<div>To print the ESTD Report, right-click in the Method Report window and select Print from the drop-down list.</div>

This page is intentionally blank.

3000 Micro GC Guide 18

Changing a Standard Module

Table 18-1 Changing a standard module

Step	Description
1	<p>A Standard GC module must be decommissioned before it can be removed. 3000 Micro GC internally communicates to each installed GC module via a unique module address.</p> <p>For a 1-2 channel or portable instrument, the module addresses are 1 and 2.</p> <p>For a 3-4 channel instrument, the module addresses are 1, 2, 3 and 4.</p> <p>Decommissioning a Standard GC module will disable the GC module’s address. Open the web browser and type the 3000 Micro GC IP Address in the address bar; press Enter to display the embedded web page.</p> <p>NOTE: At the factory, the instrument IP address is set to 10.1.1.101.</p>
2	<p>Click the Utilities tab.</p> <p>Click Change Module Configuration.</p> <div></div> <p>NOTE: Two caution messages will display. Click OK on each to continue.</p>

PN 074-558-P1A

Table 18-1 Changing a standard module (continued)

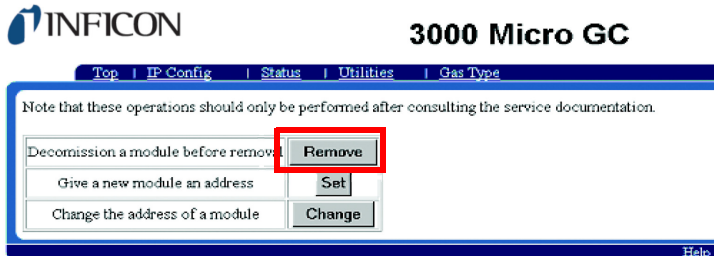
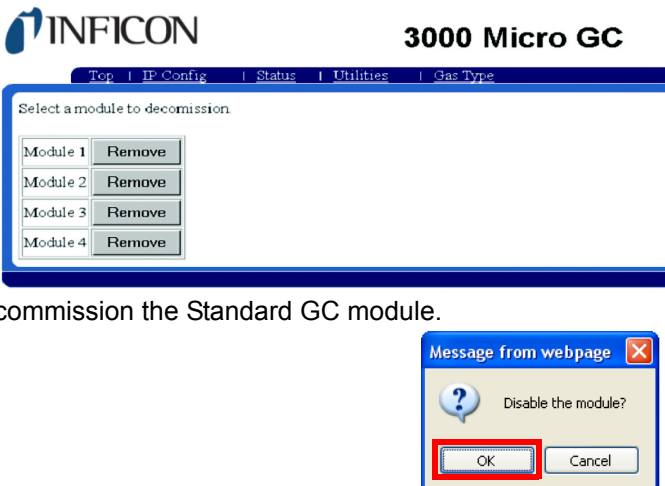





Step	Description
3	<p>Click Remove.</p> <p>NOTE: Once a Standard GC module is decommissioned, it can not be used again until it is recommissioned.</p> 
4	<p>Click Remove, adjacent to the desired Standard GC module to be decommissioned.</p>  <p>A caution will display. Click OK to decommission the Standard GC module.</p>
5	<p>Click Shutdown on the confirmation message to begin shutting down the instrument.</p> 

Table 18-1 Changing a standard module (continued)

Step	Description
6	<p>Do not turn off power. <i>Allow three minutes for the instrument to shut down.</i> A The system is shutting down... message will display.</p> <div>  <div> <p>CAUTION</p> <hr/> <p>Do not turn off the power without waiting at least three minutes. 3000 Micro GC must update configuration files before power is turned off. If power is turned off too soon, the files could be corrupted, making 3000 Micro GC unusable.</p> <hr/> </div> </div> <div>  </div>
7	<p>Set the front panel power switch to the Off position.</p>
8	<p>Remove the two screws on each side of the top cover using a Torx T-20 or Phillips Head screwdriver. Lift and remove the top cover.</p> <div>  <div> <p>WARNING - Risk Of Electric Shock</p> <hr/> <p>Electrostatic Discharge can damage electronic components. Wear a grounded wrist strap to avoid damaging 3000 Micro GC.</p> <hr/> </div> </div> <div>  </div>

PN 074-558-P1A

Table 18-1 Changing a standard module (continued)

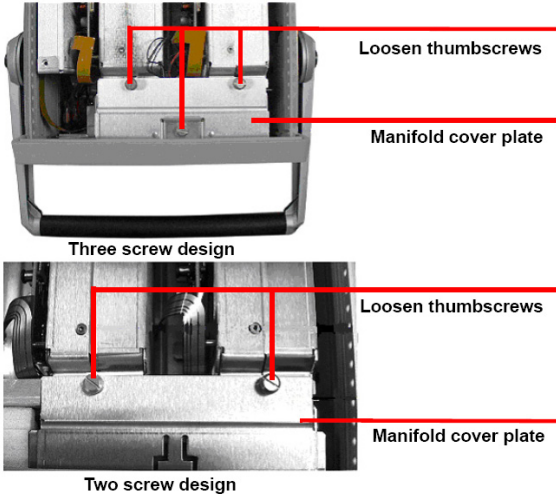
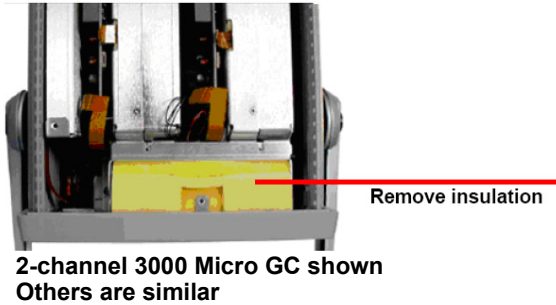
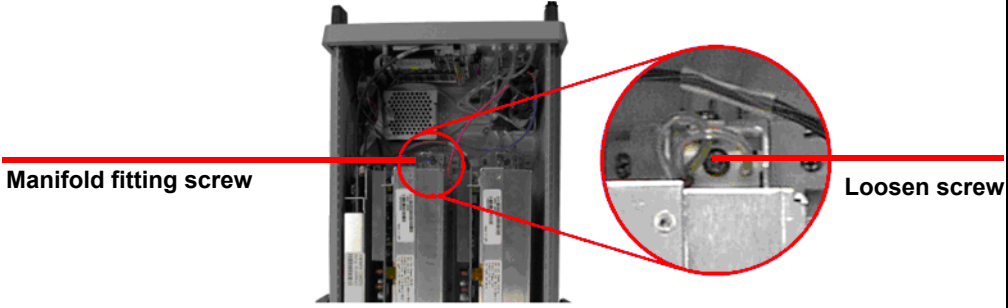

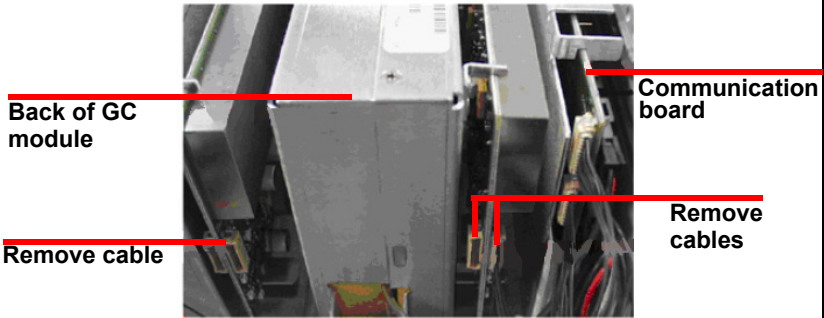
Step	Description
9	<p>Loosen the thumbscrews in the manifold cover plate.</p> <div data-bbox="820 357 1372 848">  <p>Three screw design</p> <p>Two screw design</p> </div>
10	<p>Remove the manifold cover plate. In the two screw design, slide the manifold cover plate towards the GC modules to disengage the hook in the cover plate from the tab in the chassis.</p>
11	<p>Carefully remove the manifold insulation and save it for reuse.</p> <p>NOTE: In the Portable 3000 Micro GC, the DC-DC converter assembly blocks access to the GC modules. Remove the screws that secure the DC-DC converter assembly to the side of the instrument. Without disconnecting any wires, slide the assembly off of the standoffs and gently lift it away from the GC modules.</p> <div data-bbox="820 1144 1372 1446">  <p>Remove insulation</p> <p>2-channel 3000 Micro GC shown Others are similar</p> </div>

Table 18-1 Changing a standard module (continued)

Step	Description
12	<p>Use a T-20 screwdriver (PN 5182-3468) to loosen the screw in the inlet manifold fitting at the back of the GC module.</p> <div><p>Manifold fitting screw</p><p>Loosen screw</p><p>2-channel 3000 Micro GC shown Others are similar</p></div>
13	<p>Use a 1/4 in. open-ended wrench to disconnect the inlet manifold fitting from the GC module input fitting.</p> <div><p>GC module</p><p>Inlet manifold</p><p>Disconnect inlet fittings</p><p>2-channel 3000 Micro GC shown Others are similar</p></div>
14	<p>Disconnect the communication cables leading from the GC modules (if connected) from the GC module communications board. Disconnect the power cables leading to the fan (if installed).</p> <div><p>Back of GC module</p><p>Remove cable</p><p>Communication board</p><p>Remove cables</p><p>Back view of GC modules</p></div>

PN 074-558-P1A

Table 18-1 Changing a standard module (continued)

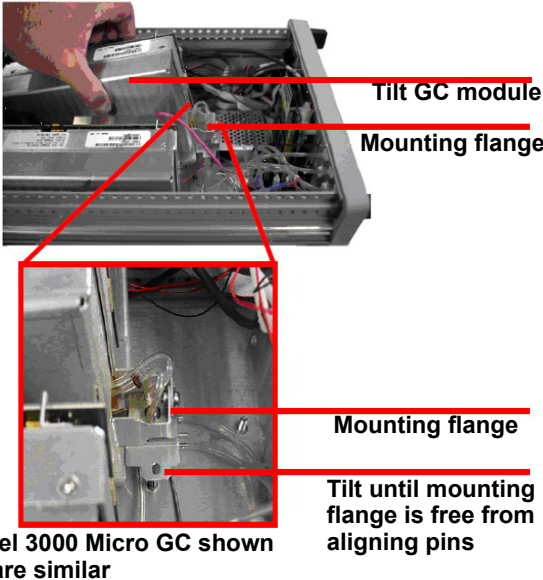

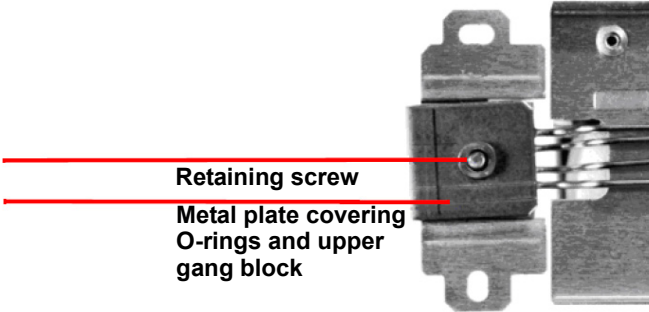
Step	Description
15	<p>Tilt the back of the GC module up until the mounting flange clears the alignment pins. Slide the GC module towards the back of the instrument until it clears the inlet manifold frame and can be lifted and removed from the front.</p>  <p>2-channel 3000 Micro GC shown Others are similar</p>
16	<p>Inspect the gang block fitting on the bottom of the chassis to make sure the mating surface is free of dust and particles.</p> 
17	<p>On the replacement module, remove the small metal plate covering the O-rings on the upper gang block.</p> 

Table 18-1 Changing a standard module (continued)

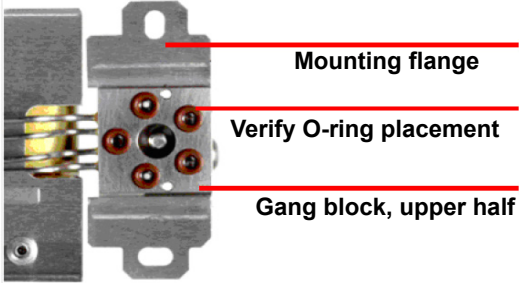
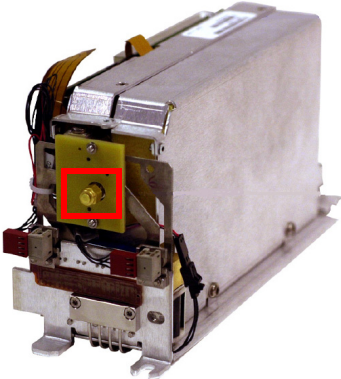
Step	Description
18	<div><div><div>Inspect the GC module mounting flange fitting to verify all new O-rings on the replacement GC module are undamaged and seated flat.</div><div><div><div>Mounting flange</div><div>Verify O-ring placement</div><div>Gang block, upper half</div></div></div></div></div>
19	<div><div><div>Use a wrench to loosen the protective 1/4 in. cap over the GC module inlet fitting.</div><div></div></div></div>

Table 18-1 Changing a standard module (continued)

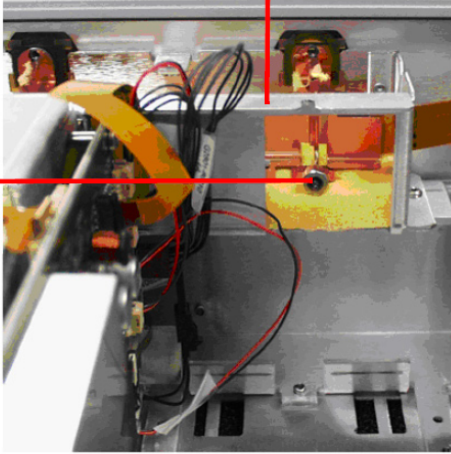

Step	Description
20	<p data-bbox="272 312 1378 417">Lower the back of the GC module into position until there is enough clearance to position it into place under the lip of the inlet manifold frame. At the same time, ensure the GC module inlet fitting mates with the Swagelok® fitting in the 3000 Micro GC inlet manifold.</p> <p data-bbox="272 428 1378 491">Slide the GC module towards the front of 3000 Micro GC until it is fully seated into the Swagelok fitting in the 3000 Micro GC inlet manifold.</p> <div data-bbox="748 533 1370 1005">  <p data-bbox="753 533 899 617">Install GC module below frame</p> <p data-bbox="753 743 899 827">Mate module manifold fittings</p> </div> <p data-bbox="997 1016 1370 1047">View from back of 3000 Micro GC</p>

Table 18-1 Changing a standard module (continued)



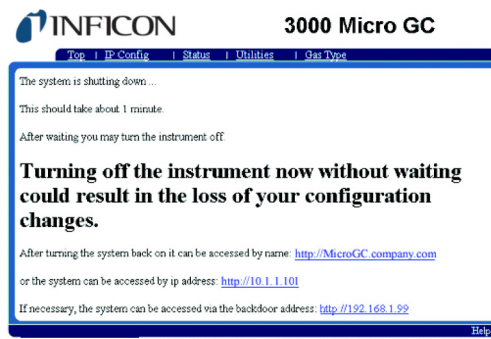
Step	Description
21	<p>Re-connect the communications cables, as shown below.</p> <p>NOTE: Do not connect more than two GC modules in series per communications board connection.</p> <p>NOTE: The GC modules and communications board use parallel communications; both connectors on each item function similarly.</p>
22	<p>Connect the inlet manifold to the GC module input fitting and tighten it finger tight. Use a 5/16 in. open-ended wrench totighten 1/4-turn past finger tight.</p>
23	<p>Tighten the screw in the mounting flange.</p>
24	<p>Carefully replace the inlet manifold insulation.</p>
25	<p>Install the manifold cover plate and the top cover.</p>
26	<p>For Standard GC modules, a new address must be given to the newly installed GC module. This can be done via the embedded web page.</p> <p>Turn the 3000 Micro GC power switch, located on its front, to the On position. The switch will illuminate green to indicate that 3000 Micro GC is on. Wait three minutes for the instrument to load.</p> <div>  <p>CAUTION</p> <p>Allow three minutes for 3000 Micro GC to update configuration files.</p> </div>

PN 074-558-P1A

Table 18-1 Changing a standard module (continued)

Step	Description
27	In the web browser address bar, type the instrument IP address.
28	Click the Utilities tab.
29	Click Change Module Configuration . <div data-bbox="812 451 1372 1018"> <p>The screenshot shows the '3000 Micro GC' web interface with the 'Utilities' tab selected. Under the 'Configuration' section, there are buttons for 'Full System', 'Module 1', 'Module 2', 'Module 3', and 'Module 4'. Below this is a 'Log' section with buttons for 'Control Module', 'Module 1', 'Module 2', 'Module 3', and 'Module 4'. At the bottom, a 'Warning!' message states: 'Before changing module configuration please consult the service documentation.' Below the warning is a button labeled 'Change Module Configuration', which is highlighted with a red box.</p> </div>
30	Click Set . <div data-bbox="560 1092 1372 1375"> <p>The screenshot shows the '3000 Micro GC' web interface with the 'Utilities' tab selected. A note states: 'Note that these operations should only be performed after consulting the service documentation.' Below the note are three buttons: 'Decommission a module before removal' (with a 'Remove' button), 'Give a new module an address' (with a 'Set' button), and 'Change the address of a module' (with a 'Change' button'). The 'Set' button is highlighted with a red box.</p> </div>
31	Use the drop-down list on the lower left portion of the screen to select the correct physical address for the new standard GC module. Each physical address corresponds to a channel number. Valid 3000 Micro GC physical addresses are 1 through 4. <div data-bbox="738 1501 1372 1774"> <p>The screenshot shows the '3000 Micro GC' web interface with the 'Utilities' tab selected. Under the 'Adding a new module' section, there is a 'Select a new address:' label and a note: 'Note that only 1 and 2 are currently supported addresses'. Below this is a drop-down list with the following options: 1, 2, 3, 4, 5, 6, 7, 8. The value '1' is selected and highlighted with a red box.</p> </div>

Table 18-1 Changing a standard module (continued)



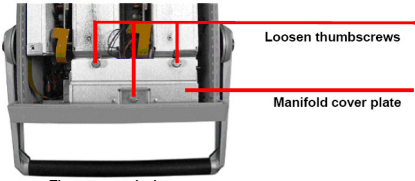
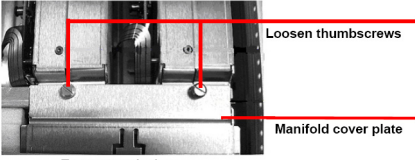
Step	Description
32	<p>After selecting the address, click Restart.</p> 
33	<p>A The system is shutting down... message is displayed. Allow three minutes for the instrument to update configuration files.</p> <div data-bbox="373 724 446 787">  </div> <div data-bbox="467 735 625 772"> <h3>CAUTION</h3> </div> <div data-bbox="490 808 1295 934"> <p>Do not turn off the power without waiting at least three minutes. 3000 Micro GC must update configuration files before power is turned off. If power is turned off too soon, the files could be corrupted and render 3000 Micro GC unusable.</p> </div> 
34	<p>Open the web browser and type the 3000 Micro GC IP Address in the address bar. Verify that the recently installed replacement module is available in the module list.</p>
35	<p>Once the replacement standard module is installed, update EZ IQ with the revised 3000 Micro GC hardware configuration. For detailed instructions on how to configure EZ IQ, see 074-537 EZ IQ Installation Guide.</p>

This page is intentionally blank.

3000 Micro GC Guide 19

Changing a Performance Enhanced Module

Table 19-1 Changing a Performance Enhanced module

Step	Description
1	<p>To replace a Performance Enhanced module with another Performance Enhanced module, turn of the instrument by setting the front panel power switch to the Off position. Remove the two screws on each side of the top cover using a Torx T-20 or Phillips Head screwdriver. Lift and remove the top cover.</p> <div><div></div><div>WARNING - Risk Of Electric Shock Electrostatic Discharge can damage electronic components. Wear a grounded wrist strap to avoid damaging 3000 Micro GC.</div></div> <div></div>
2	<p>Loosen the thumbscrews in the manifold cover plate.</p> <div><div> Three screw design</div><div> Two screw design</div></div>
3	<p>Remove the manifold cover plate. Slide the manifold cover plate towards the GC module to disengage the hook in the cover plate from the tab in the chassis.</p>

PN 074-558-P1A

Table 19-1 Changing a Performance Enhanced module (continued)

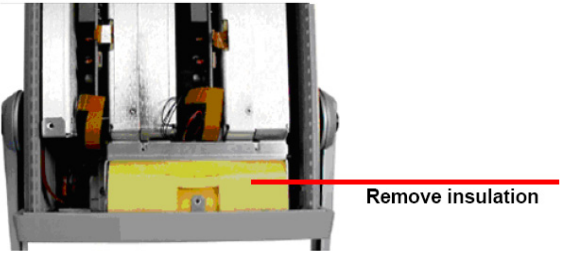
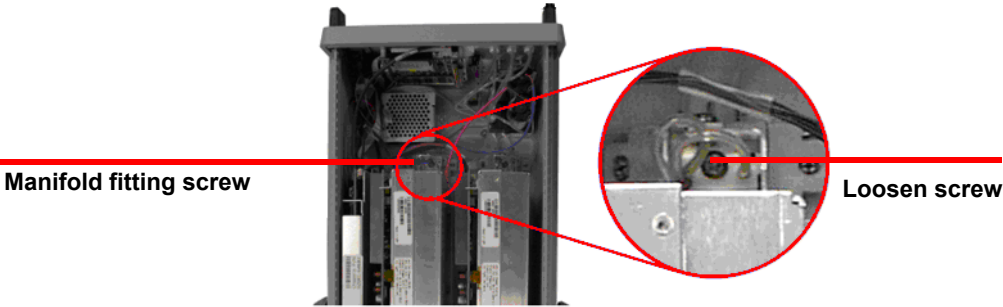
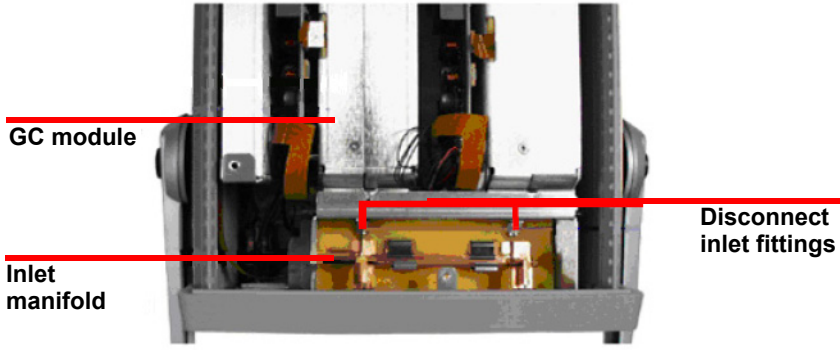
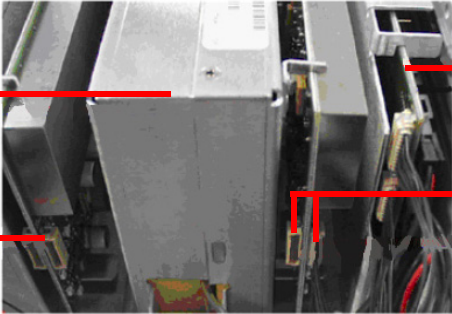
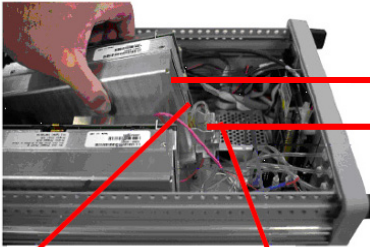
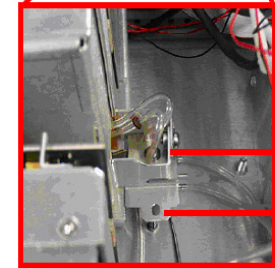
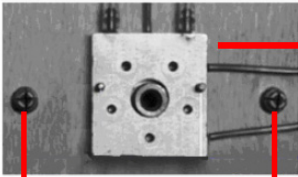
Step	Description
4	<p>Carefully remove the manifold insulation. Save for reuse.</p> <p>NOTE: In the Portable 3000 Micro GC, the DC-DC converter assembly blocks access to the GC modules. Remove the screws that secure the DC-DC converter assembly to the side of the instrument. Without disconnecting any wires, slide the assembly off of the standoffs and gently lift it away from the GC modules.</p>  <p>Remove insulation</p> <p>2-channel 3000 Micro GC shown Others are similar</p>
5	<p>Use a T-20 screwdriver to loosen the screw in the inlet manifold fitting at the back of the GC module.</p>  <p>Manifold fitting screw</p> <p>Loosen screw</p> <p>2-channel 3000 Micro GC shown</p>
6	<p>Use a 1/4 in. open-ended wrench to disconnect the inlet manifold fitting from the GC module input fitting.</p>  <p>GC module</p> <p>Inlet manifold</p> <p>Disconnect inlet fittings</p> <p>2-channel 3000 Micro GC shown Others are similar</p>

Table 19-1 Changing a Performance Enhanced module (continued)

Step	Description
7	<p>Disconnect the communication cables leading from the GC modules, if present, to the communications board connectors for the GC modules. If needed, also disconnect the power cables leading to the fan.</p> <div><div><div>Back of GC module</div><div>Remove cable</div></div><div><div>Communication board</div><div>Remove cables</div></div><div>Back view of GC modules</div></div>
8	<div><div><div><div><div>Tilt GC module</div></div><div><div>Mounting flange</div></div></div><div>2-channel 3000 Micro GC shown Others are similar</div><div>Tilt until mounting flange is free from aligning pins</div></div><p>Tilt the back of the GC module up until the mounting flange clears the alignment pins. Slide the GC module towards the back of the instrument until it clears the inlet manifold frame and can be lifted and removed from the front.</p></div>
9	<p>Inspect the gang block fitting on the bottom of the chassis to ensure the mating surface is clean.</p> <div><div><div>Gang block, lower half</div></div><div>Alignment pins</div></div>

PN 074-558-P1A

Table 19-1 Changing a Performance Enhanced module (continued)

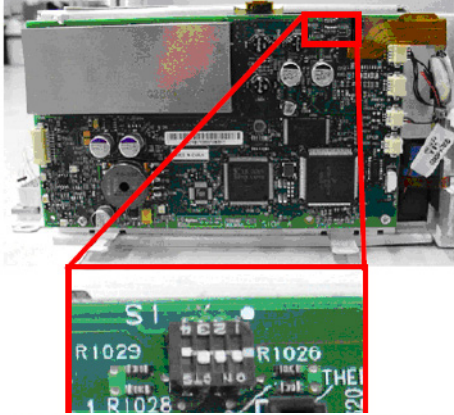

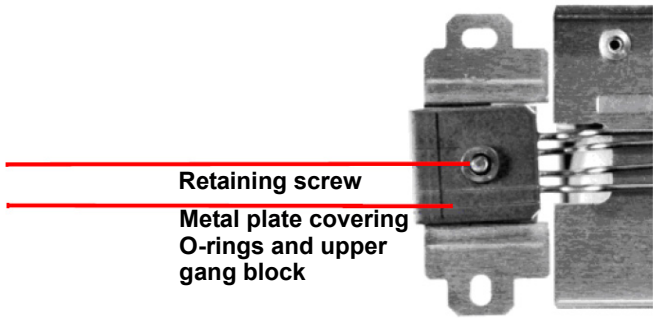
Step	Description
10	<p data-bbox="272 310 1377 380">On the replacement module, set the Performance Enhanced GC module address by setting the DIP switch on the upper right of the D-board to the appropriate configuration.</p> <p data-bbox="272 394 1377 464">Each setting determines a unique module address. Please refer to the label on the top of the module. The DEF setting is the factory default DIP switch setting.</p> <div data-bbox="857 478 1308 890">  </div> <div data-bbox="800 894 1382 1245">  </div>
11	<p data-bbox="272 1266 1170 1299">Remove the small metal plate covering the O-rings on the upper gang block.</p> <div data-bbox="703 1329 1352 1644">  <p data-bbox="878 1493 1062 1522">Retaining screw</p> <p data-bbox="878 1530 1109 1612">Metal plate covering O-rings and upper gang block</p> </div>

Table 19-1 Changing a Performance Enhanced module (continued)

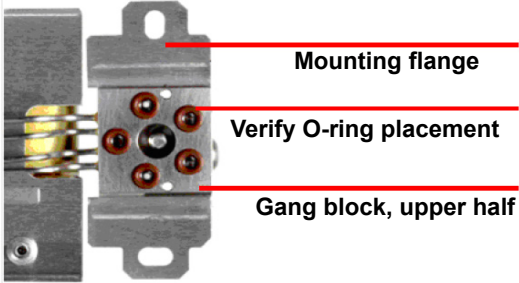
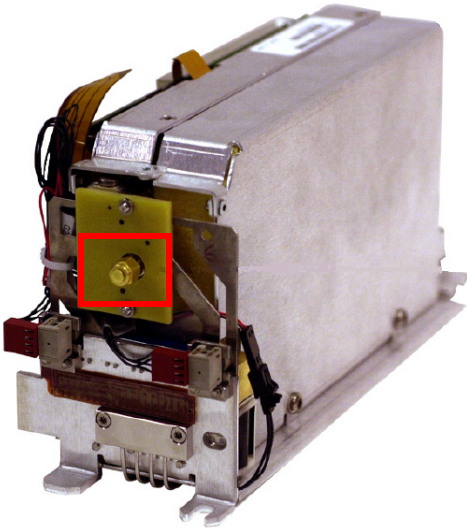
Step	Description
12	<div><div><div>Inspect the GC module mounting flange fitting to verify all new O-rings on the replacement GC module are undamaged and seated flat.</div><div><div><div>Mounting flange</div><div>Verify O-ring placement</div><div>Gang block, upper half</div></div></div></div></div>
13	<div><div><div>Remove the protective 1/4 in. cap over the GC module inlet fitting.</div><div></div></div></div>

Table 19-1 Changing a Performance Enhanced module (continued)

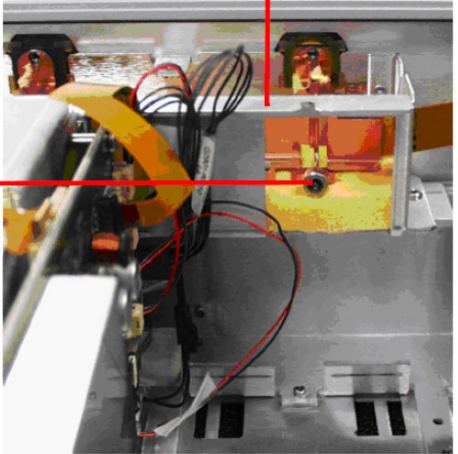
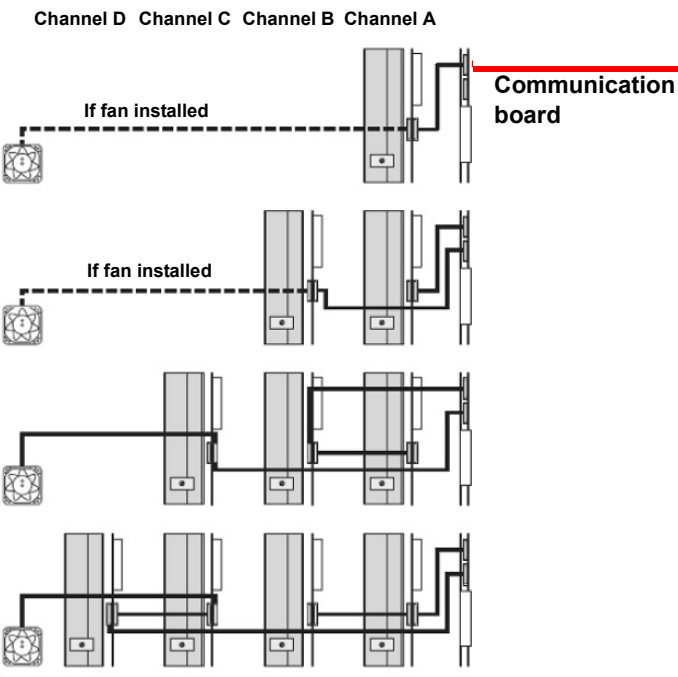
Step	Description
14	<p data-bbox="272 310 1377 415">Lower the back of the GC module into position until there is enough clearance to position it into place under the lip of the inlet manifold frame. At the same time, ensure the GC module inlet fitting mates with the Swagelok fitting in the 3000 Micro GC inlet manifold.</p> <p data-bbox="272 426 1377 499">Slide the GC module towards the front of 3000 Micro GC until it is fully seated into the Swagelok fitting in the 3000 Micro GC inlet manifold.</p> <div data-bbox="751 531 1377 1003"> <p data-bbox="751 531 898 615">Install GC module below frame</p> <p data-bbox="751 741 898 825">Mate module manifold fittings</p>  </div> <p data-bbox="995 1014 1377 1045">View from back of 3000 Micro GC</p>

Table 19-1 Changing a Performance Enhanced module (continued)

Step	Description
15	<p>Re-connect the communications cables.</p> <p>NOTE: Do not connect more than two GC modules in series per communications board connection.</p> <p>NOTE: The GC modules and communications board use parallel communications; both connectors on each item function similarly.</p> 
16	Connect the inlet manifold to the GC module input fitting and tighten it finger tight. Use a 5/16 in. open-ended wrench, to tighten 1/4-turn past finger tight.
17	Tighten the screw in the mounting flange.
18	Carefully replace the inlet manifold insulation.
19	Install the manifold cover plate and the top cover.
20	<p>Turn the 3000 Micro GC power switch, located on its front, to the on position. The switch will illuminate green to indicate that 3000 Micro GC is on.</p> <p>Wait three minutes for the instrument to load.</p>
21	Open the web browser and type the 3000 Micro GC IP Address in the address bar. Verify that the newly installed replacement module is available in the module list.
22	Once the replacement GC module is installed, update EZ IQ with the updated 3000 Micro GC hardware configuration. For detailed instructions on how to configure EZ IQ, consult the PN 074-537-P1A EZ IQ Installation Guide.

This page is intentionally blank.

3000 Micro GC Guide 20

Replacing a Standard Module with a Performance Enhanced Module

Table 20-1 Changing a standard module

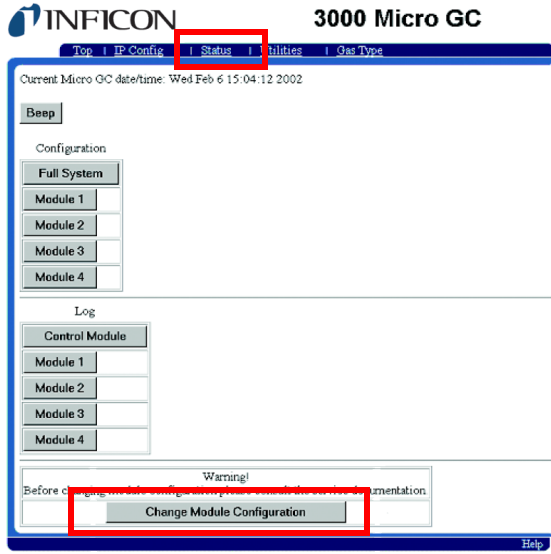
Step	Description
1	<p>A Standard GC module must be decommissioned before it can be removed. 3000 Micro GC internally communicates to each installed GC module using a unique module address.</p> <p>For a 1-2 channel or portable instrument, the module addresses are 1 and 2.</p> <p>For a 3-4 channel instrument, the module addresses are 1, 2, 3 and 4.</p> <p>Decommissioning a Standard GC module will disable the GC module's address.</p> <p>Open the web browser and type the 3000 Micro GC IP Address in the address bar, then press Enter to display the embedded web page.</p> <p>NOTE: At the factory, the instrument IP address is set to 10.1.1.101.</p>
2	<p>Click the Utilities tab.</p> <p>Click Change Module Configuration.</p> 
3	<p>Two caution messages will display. Click OK on each to continue.</p>

Table 20-1 Changing a standard module (continued)

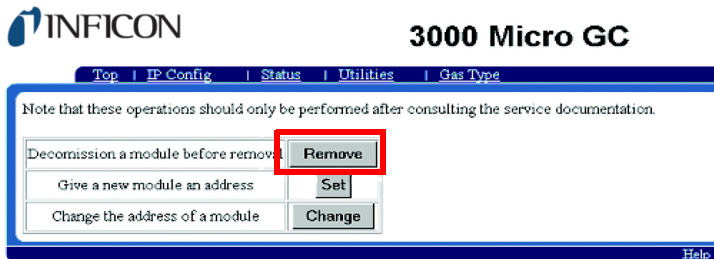




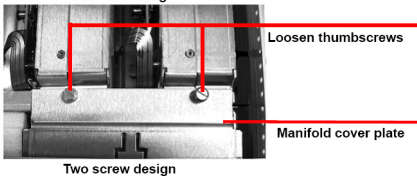
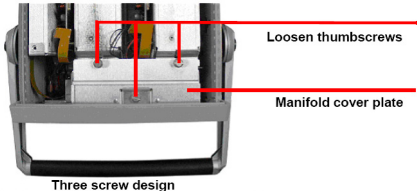
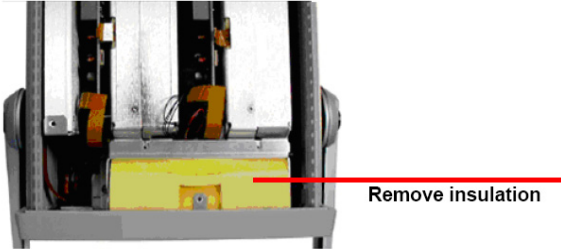
Step	Description
4	<p>Click Remove.</p> <p>NOTE: Once a Standard GC module is decommissioned, it cannot be used until it is recommissioned.</p>  <p>The screenshot shows the INFICON 3000 Micro GC interface. At the top, there is a navigation bar with links: Top, IP Config, Status, Utilities, and Gas Type. Below the navigation bar, a note states: 'Note that these operations should only be performed after consulting the service documentation.' Underneath the note, there are three buttons: 'Remove', 'Set', and 'Change'. The 'Remove' button is highlighted with a red box.</p>
5	<p>Click Remove next to the Standard GC module to decommission. A caution will display. Click OK to decommission the Standard GC module.</p>  <p>The screenshot shows the INFICON 3000 Micro GC interface. At the top, there is a navigation bar with links: Top, IP Config, Status, Utilities, and Gas Type. Below the navigation bar, the text 'Select a module to decommission' is displayed. Underneath, there are four rows, each with a module number and a 'Remove' button. The 'Remove' button for Module 1 is highlighted with a red box.</p>
6	<p>Click Shutdown on the confirmation message. Allow three minutes for the instrument to shut down.</p>  <p>The screenshot shows the INFICON 3000 Micro GC interface. At the top, there is a navigation bar with links: Top, IP Config, Status, Utilities, and Gas Type. Below the navigation bar, the text 'Module 1 has been decommissioned' is displayed. Underneath, there is a 'Shutdown' button, which is highlighted with a red box.</p>
7	<p>Set the front panel power switch to the Off position.</p>

Table 20-1 Changing a standard module (continued)

Step	Description
8	<p>Remove the two screws on each side of the top cover using a Torx T-20 or Phillips Head screwdriver. Lift and remove the top cover.</p> <div>WARNING - Risk Of Electric Shock<p>Electrostatic Discharge can damage electronic components. Wear a grounded wrist strap to avoid damaging 3000 Micro GC.</p></div> <div></div>
9	<p>Loosen the thumbscrews in the manifold cover plate.</p> <div></div>
10	<p>Remove the manifold cover plate. In the two screw design, slide the manifold cover plate towards the GC module to disengage the hook in the cover plate from the tab in the chassis.</p>
11	<p>Carefully remove the manifold insulation. Save it for reuse.</p> <p>NOTE: In the Portable 3000 Micro GC, the DC-DC converter assembly blocks access to the GC modules. Remove the screws that secure the DC-DC converter assembly to the side of the instrument. Without disconnecting any wires, slide the assembly off of the standoffs and gently lift it away from the GC modules.</p> <div><p>2-channel 3000 Micro GC shown Others are similar</p></div>

PN 074-558-P1A

Table 20-1 Changing a standard module (continued)

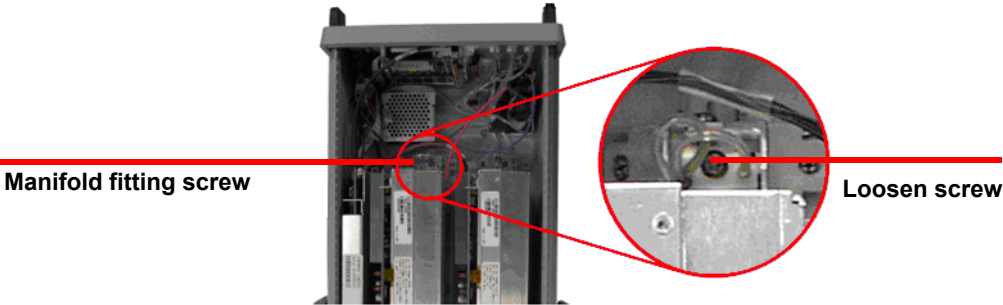
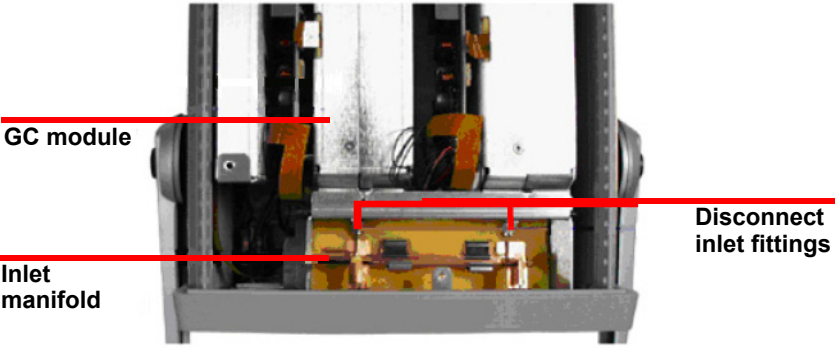
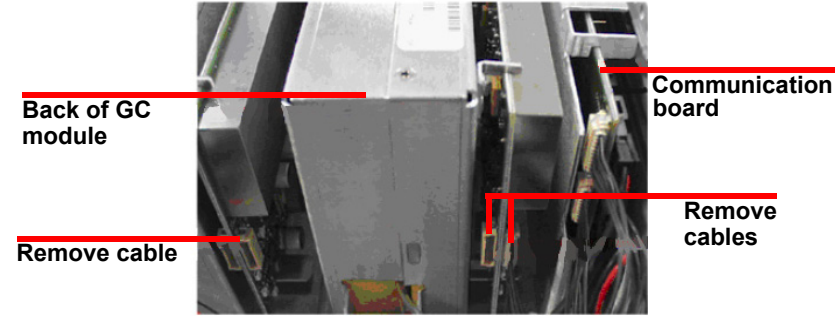
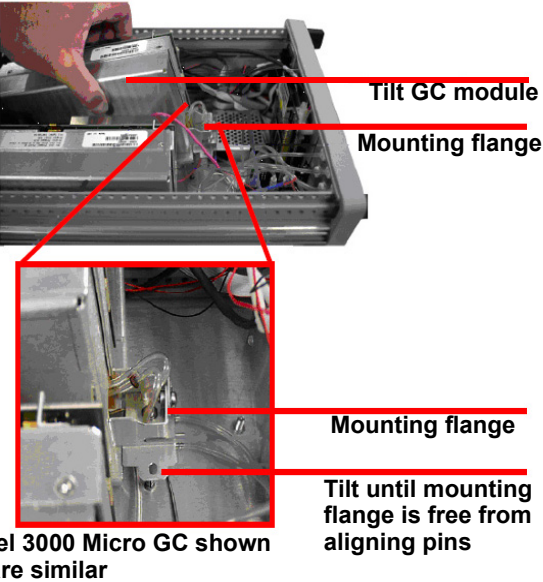
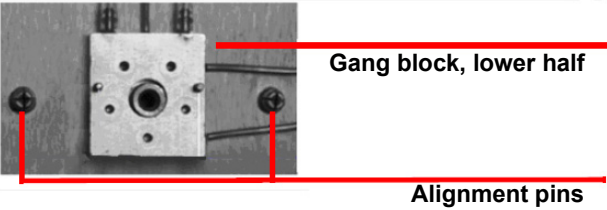
Step	Description
12	<p>Use a T-20 screwdriver to loosen the screw in the inlet manifold fitting at the back of the GC module.</p>  <p>Manifold fitting screw</p> <p>Loosen screw</p> <p>2-channel 3000 Micro GC shown</p>
13	<p>Use a 1/4 in. open-ended wrench to disconnect the inlet manifold fitting from the GC module input fitting.</p>  <p>GC module</p> <p>Inlet manifold</p> <p>Disconnect inlet fittings</p> <p>2-channel 3000 Micro GC shown Others are similar</p>
14	<p>Disconnect the communication cables leading from the GC modules, if present, to the communications board connectors for the GC modules. If needed, also disconnect the power cables leading to the fan.</p>  <p>Back of GC module</p> <p>Remove cable</p> <p>Communication board</p> <p>Remove cables</p> <p>Back view of GC modules</p>

Table 20-1 Changing a standard module (continued)

Step	Description
15	<p>Tilt the back of the GC module up until the mounting flange clears the alignment pins. Slide the GC module towards the back of the instrument until it clears the inlet manifold frame and can be lifted and removed from the front.</p> <div><p>Tilt GC module</p><p>Mounting flange</p><p>Mounting flange</p><p>Tilt until mounting flange is free from aligning pins</p><p>2-channel 3000 Micro GC shown Others are similar</p></div>
16	<p>Inspect the gang block fitting on the bottom of the chassis to make sure the mating surface is free of dust and particles.</p> <div><p>Gang block, lower half</p><p>Alignment pins</p></div>

PN 074-558-P1A

Table 20-1 Changing a standard module (continued)

Step	Description
17	<p>On the new module, set the Performance Enhanced GC module address by setting the DIP switch on the upper right of the D-board to the appropriate configuration. Each setting determines a unique module address. Please refer to the label on the top of the module. The DEF setting is the factory default DIP switch setting.</p> <div data-bbox="857 457 1308 894" data-label="Image"> </div> <div data-bbox="800 919 1382 1272" data-label="Image"> </div>
18	<p>Remove the small metal plate covering the O-rings on the upper gang block.</p> <div data-bbox="704 1356 1352 1671" data-label="Image"> </div>

Table 20-1 Changing a standard module (continued)

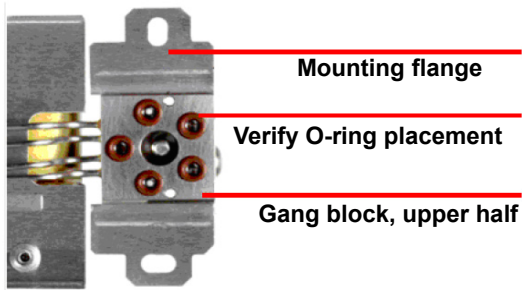
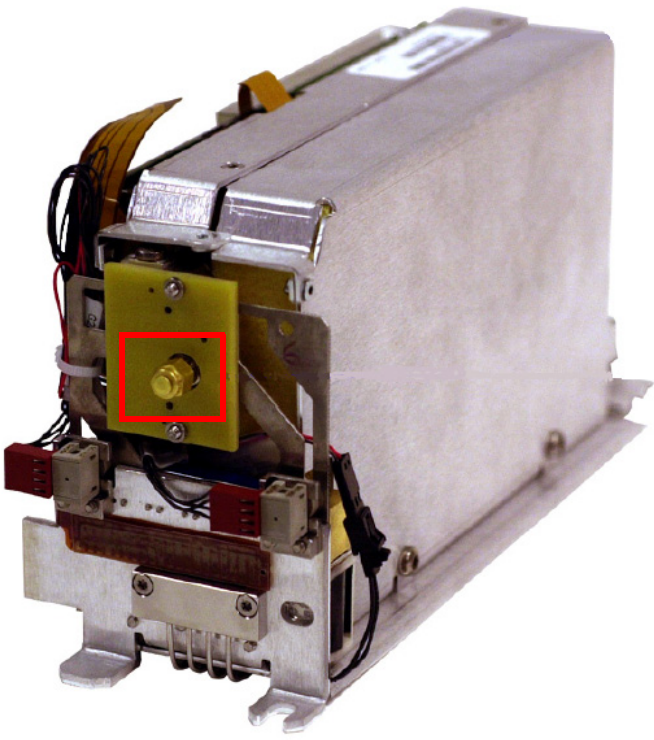
Step	Description
19	<p>Inspect the GC module mounting flange fitting to verify all new O-rings on the replacement GC module are undamaged and seated flat.</p> <div></div>
20	<p>Remove the protective 1/4 in. cap over the GC module inlet fitting.</p> <div></div>

Table 20-1 Changing a standard module (continued)

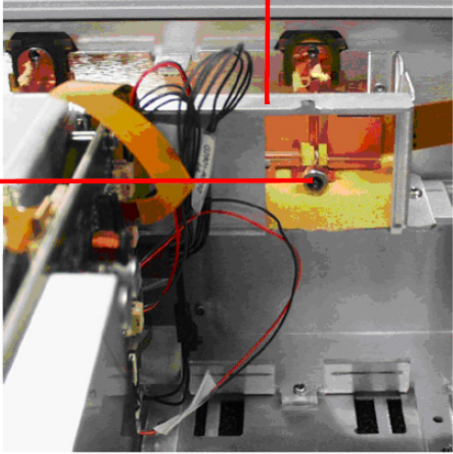
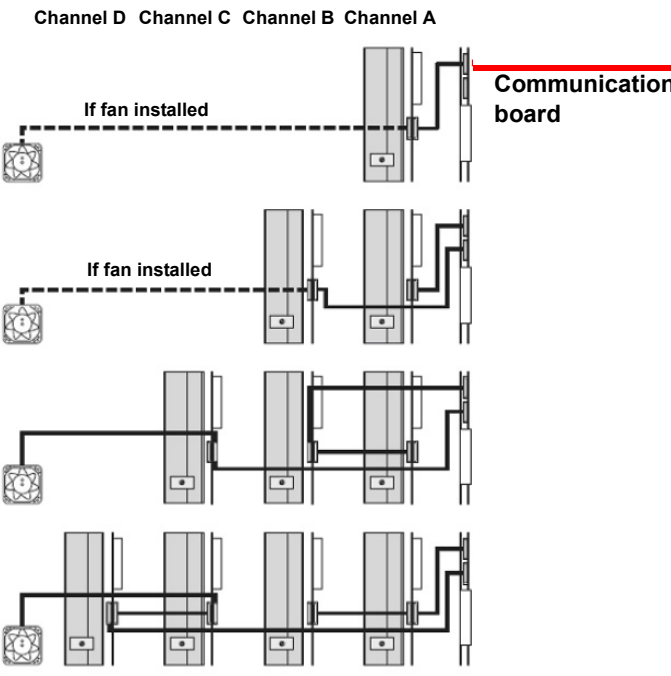
Step	Description
21	<p data-bbox="272 312 1372 417">Lower the back of the GC module into position until there is enough clearance to position it into place under the lip of the inlet manifold frame. At the same time, ensure the GC module inlet fitting mates with the Swagelok fitting in the 3000 Micro GC inlet manifold.</p> <p data-bbox="272 428 1372 491">Slide the GC module towards the front of the 3000 Micro GC until it is fully seated into the Swagelok fitting in the 3000 Micro GC inlet manifold.</p> <div data-bbox="748 533 1372 1005">  <p data-bbox="753 533 894 617">Install GC module below frame</p> <p data-bbox="753 743 902 827">Mate module manifold fittings</p> </div> <p data-bbox="997 1016 1372 1047">View from back of 3000 Micro GC</p>

Table 20-1 Changing a standard module (continued)

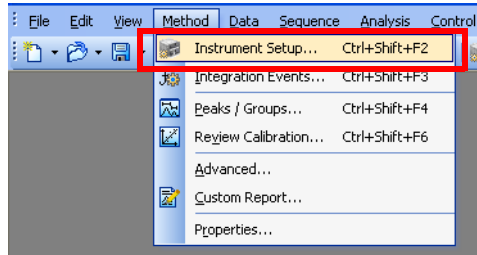
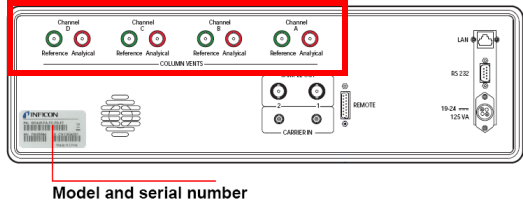
Step	Description
22	<p>Re-connect the communications cables.</p> <p>NOTE: Do not connect more than two GC modules in series per communications board connection.</p> <p>NOTE: The GC modules and communications board use parallel communications; both connectors on each item function similarly.</p> 
23	Connect the inlet manifold to the GC module input fitting and tighten it finger tight. Using a 5/16 in. open-ended wrench, tighten 1/4-turn past finger tight.
24	Tighten the screw in the mounting flange.
25	Carefully replace the inlet manifold insulation.
26	Install the manifold cover plate and the top cover.
27	Turn the 3000 Micro GC power switch, located on its front, to the on position. The switch will illuminate green to indicate that 3000 Micro GC is on. Wait three minutes for the instrument to load.
28	Open the web browser and type the 3000 Micro GC IP Address in the address bar. Verify that the recently installed replacement module is available in the module list.
29	After installing the replacement GC module, update EZ IQ with the updated 3000 Micro GC hardware configuration. For detailed instructions on how to configure EZ IQ, consult the PN 074-537 EZ IQ Installation Guide.

This page is intentionally blank.

3000 Micro GC Guide 21

Testing Instrument Flow Rates

Table 21-1 Test instrument flow rates

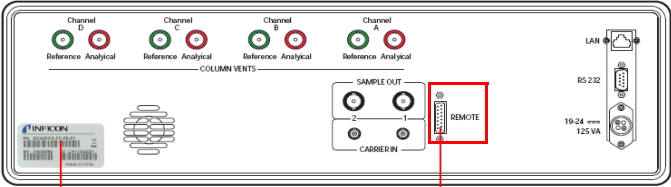
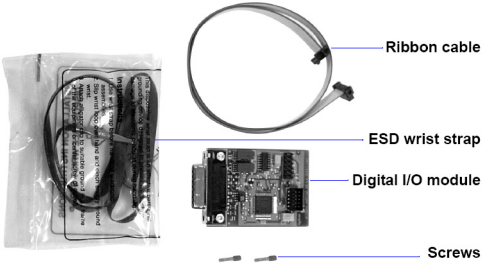
Step	Description
1	<p>Instrument flow rates can be tested for diagnostic purposes. Prior to testing, create a test method in EZ IQ.</p> <p>Click File >> Method >> Create.</p>
2	<p>Click Method >> Instrument Setup.</p> 
3	<p>In the Instrument Setup window, set the Column Head Pressure and Post Column Pressure to 20 PSI.</p>
4	<p>Download the method by clicking Control >> Download.</p>
5	<p>Using a flowmeter, measure the flow rates from the Analytical and Reference column vents, located on the back of 3000 Micro GC. Male luer locks are provided in the instrument ship kit. The female end will attach directly onto the column vents. The ship kit also contains a 1/16 in. piece of polyurethane tubing, which can be attached to the male end.</p> <p>Analytical flow rate should be 1.0 +/- 0.5 mL/min.</p> <p>Reference flow rate should be 0.6 +/- 0.4 mL/min.</p> <p>NOTE: Flow rates will vary depending on column length, internal diameter, and film thickness.</p> 
6	<p>If no flowmeter is available, a rudimentary test can be performed to verify the presence of flow:</p> <ol style="list-style-type: none"> 1 Attach a male luer lock to the piece of 1/16 inch polyurethane tubing, provided in the instrument shipping kit. 2 Attach the female end of the luer lock to one of the column vents on the back of the instrument. Set the open end of the tubing in a container of water. Bubbles indicate flow.
3	<p>If there is no flow from either or both the Analytical and Reference ports, contact INFICON.</p>

This page is intentionally blank.

3000 Micro GC Guide 22

Installing the Digital I/O Module

Table 22-1 Installing the digital I/O module

Step	Description
1	<p>For some applications, an external triggering device may be required for sampling. Install the Digital I/O Module (PN G2847A) in 3000 Micro GC and connect it to the external triggering device via a Remote Start Cable (PN G2801-60618).</p> <p>NOTE: A 4-channel 3000 Micro GC is displayed below.</p> <div><p>Model and serial number</p><p>Input to Digital I/O Module</p></div>
2	<p>The digital I/O kit contains the following items:</p> <div><p>Ribbon cable</p><p>ESD wrist strap</p><p>Digital I/O module</p><p>Screws</p></div>
3	<p>The following tools are required to install the digital I/O module (not provided):</p> <ul style="list-style-type: none">• Torx T-10 driver• Torx T-20 screwdriver driver• Needle nose pliers (helpful for disconnecting cables)• Phillips Head Screwdriver• Flat-bladed screwdriver

PN 074-558-P1A

Table 22-1 Installing the digital I/O module (continued)

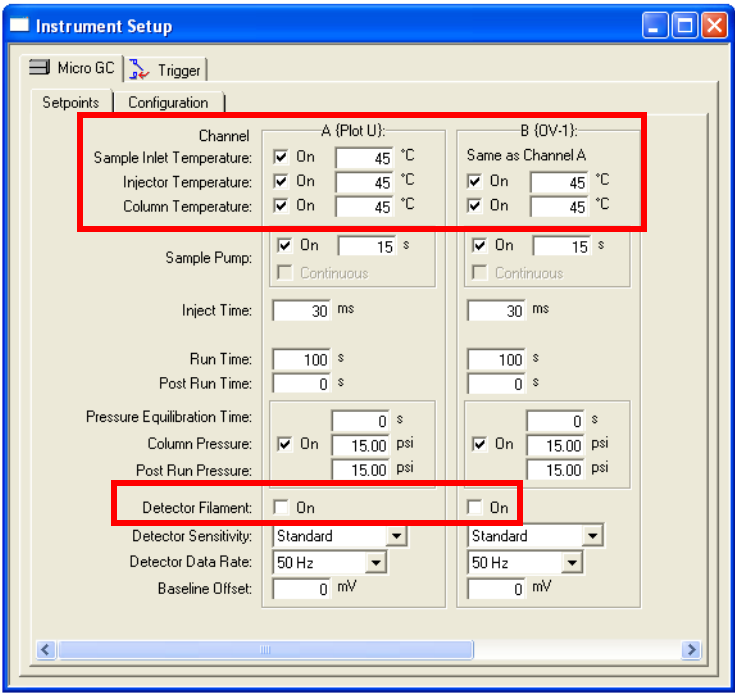




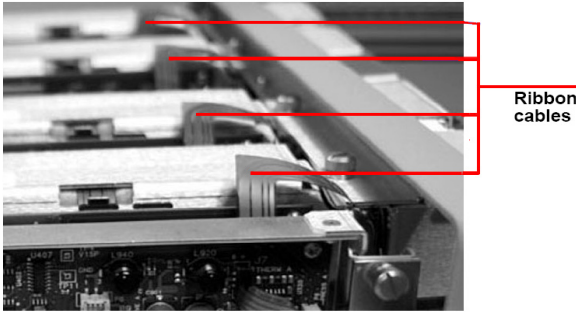
Step	Description
4	<p>Before installing the digital I/O module, prepare the instrument for service. Lower the heated inlets, injectors, and columns to 45°C. Clear Detector Filament: to turn off the detectors.</p> 
5	<p>Download the method from EZ IQ onto 3000 Micro GC. Check the EZ IQ Instrument Status screen by clicking Control >> Instrument Status. Wait until the method setpoints are reached (i.e., the instrument is in the Ready state).</p>
6	<p>Exit EZ IQ.</p>
7	<p>Set the front panel power switch to the off position. Disconnect the 3000 Micro GC power supply from the back of the instrument.</p> <div>  <div> <p>WARNING</p> <hr/> <p>During this process, the 3000 Micro GC internal components will be exposed. To avoid damaging 3000 Micro GC, turn the power switch off and disconnect all external power to 3000 Micro GC.</p> <hr/> </div> </div>

Table 22-1 Installing the digital I/O module (continued)

Step	Description
8	<p>Remove the two screws on each side of the top cover using a Torx T-20 or Phillips Head screwdriver. Lift and remove the top cover.</p> <div><div></div><div>WARNING - Risk Of Electric Shock Electrostatic Discharge can damage electronic components. Wear a grounded wrist strap to avoid damaging 3000 Micro GC.</div></div> <div><div>Screws</div></div>
9	<p>Unscrew and remove the four black feet from the back panel using a Torx 20 screwdriver. Retain the screws for reassembly. Carefully slide the main chassis out of the back of the 3000 Micro GC until the communication board is exposed.</p> <div><div></div><div>CAUTION Do not damage the ribbon cable near the upper side portion of each module.</div></div> <div></div>
10	<p>If removing a digital I/O module, refer to step 11.</p> <p>If installing a digital I/O module, follow steps 13 through 23.</p>

PN 074-558-P1A

Table 22-1 Installing the digital I/O module (continued)

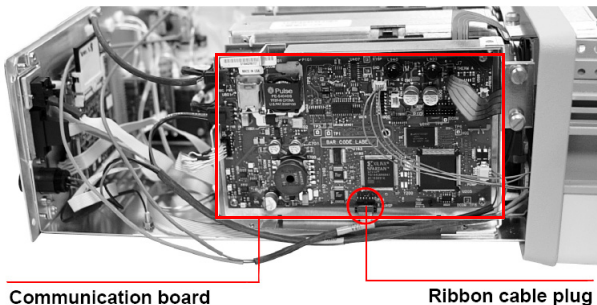
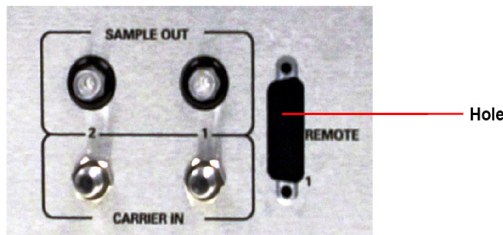
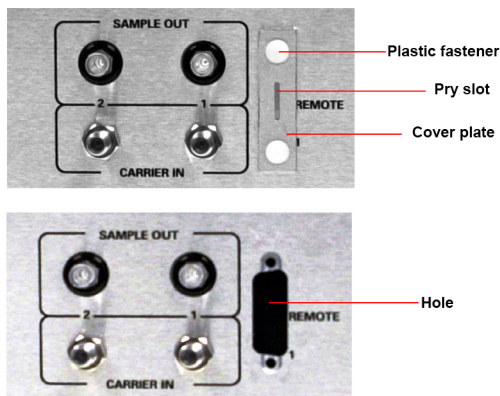
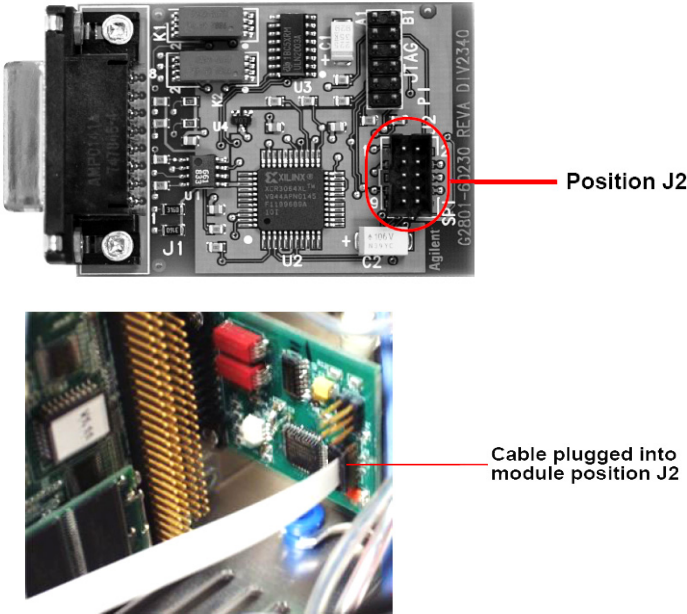
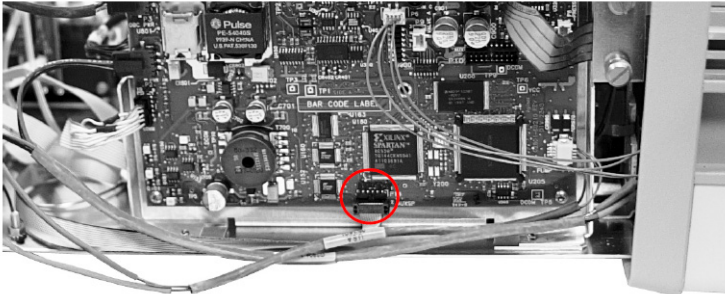
Step	Description
11	<p>Disconnect the ribbon cable from the communication board.</p> 
12	<p>Remove the two screws fastening the module to the back panel. Retain the screws for later use. Remove the digital I/O module and ribbon cable.</p> 
13	<p>To install a digital I/O module, use a flat-head screwdriver to pry the cover plate off the instrument's back panel.</p> 
14	<p>Slide the new Digital I/O Module through the hole in the back panel so that the two threaded studs protrude through. Use the retained screws to secure it in place.</p>

Table 22-1 Installing the digital I/O module (continued)




Step	Description
15	<p>Attach one end of the ribbon cable to digital I/O module position J2.</p> 
16	<p>Attach the other end of the cable to the communication board at position P14.</p> 
17	<p>Carefully slide the chassis back into the instrument. Take care not to damage any wiring or cables. Use the retained screws to reassemble the top cover. Attach all feet.</p>
18	<p>Turn the 3000 Micro GC power switch, located on its front, to the On position. The switch will illuminate green to indicate that 3000 Micro GC is on.</p> <p>Wait three minutes for the instrument to load.</p>
19	<p>Refer to the 074-519 3000 Micro GC Operating Manual for details on how to use the digital I/O module.</p>

PN 074-558-P1A

This page is intentionally blank.

3000 Micro GC Guide 23 Replacing the Battery Cable Fuse on a Portable Instrument

Table 23-1 Replacing the battery cable fuse on a portable instrument

Step	Description
1	Before beginning work on an instrument, use EZ IQ software to prepare the instrument for service. Lower the heated inlets, injectors, and columns to 45°C. Turn off the Detector Filaments .
2	Download the method from EZ IQ onto 3000 Micro GC. Check the EZ IQ Instrument Status screen by clicking Control >> Instrument Status . Wait until the method setpoints are reached (i.e., the instrument is in the Ready state).
3	Shut down EZ IQ.
4	<p>Set the front panel power switch to the Off position. Disconnect the 3000 Micro GC power supply from the back of the instrument.</p> <div>  <p>WARNING</p> <p>During this process, the 3000 Micro GC internal components will be exposed. To avoid damaging 3000 Micro GC, turn the power switch off and disconnect all external power to 3000 Micro GC.</p> </div>
5	<p>Remove the two screws on each side of the top cover, using a Torx T-20 screwdriver or a Phillips Head screwdriver. Lift and remove the top cover.</p> <div>  <p>WARNING - Risk Of Electric Shock</p> <p>Electrostatic Discharge can damage electronic components. Wear a grounded wrist strap to avoid damaging 3000 Micro GC.</p> </div> <div>  </div>

PN 074-558-P1A

Table 23-1 Replacing the battery cable fuse on a portable instrument (continued)

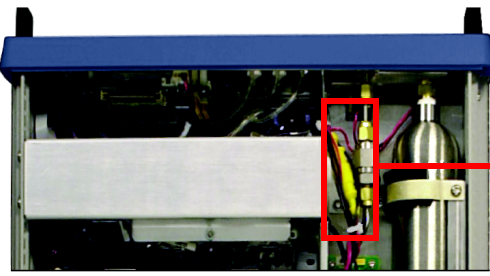
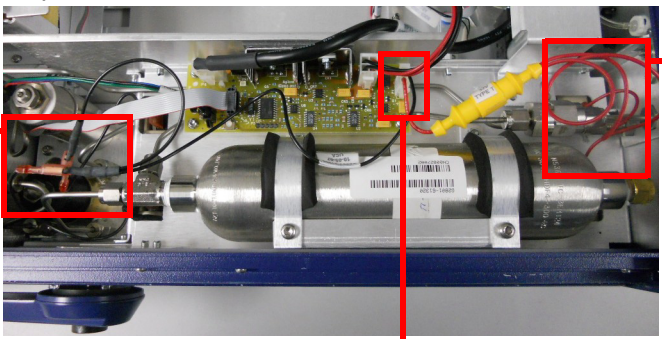
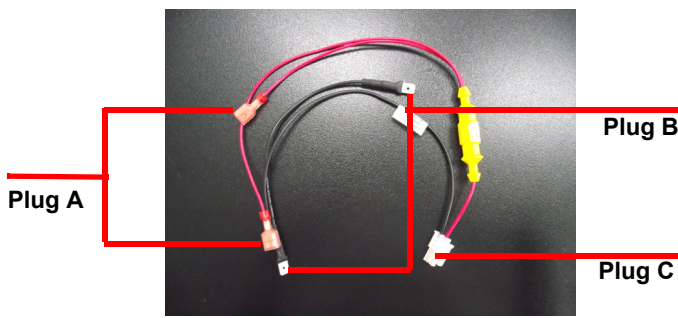
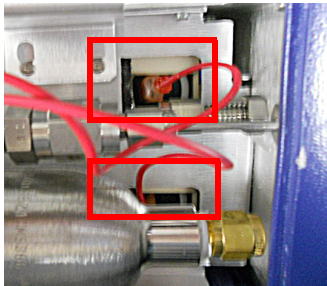
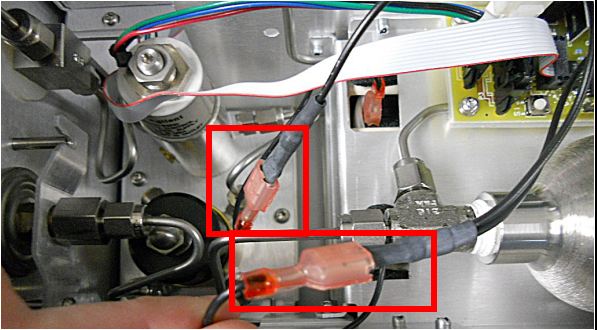
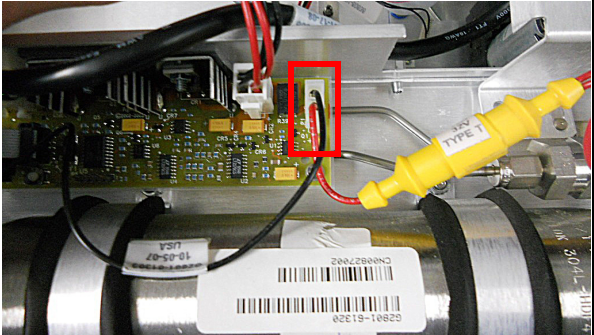
Step	Description
6	<p>Locate the battery cable fuse.</p>  <p>Fuse, 3AG Slow-blow, 10 A 32 V Type T</p>
7	<p>Locate the battery cable fuse positions.</p>  <p>Position 1</p> <p>Position 2</p> <p>Position 3</p>
8	<p>Disconnect the battery cable fuse connectors in Position 1, Position 2, and Position 3.</p>
9	<p>Remove the used battery cable fuse. Locate Plug A, Plug B, and Plug C.</p> <p>NOTE: The battery cable fuse contains two Plug A connectors, two Plug B connectors, and one Plug C connector.</p>  <p>Plug A</p> <p>Plug B</p> <p>Plug C</p>
10	<p>On the replacement battery cable fuse, connect both Plug A connectors to Position 1.</p> 

Table 23-1 Replacing the battery cable fuse on a portable instrument (continued)

Step	Description
11	<div>Next, connect both Plug B connectors to Position 3.</div> <div></div>
12	<div>Connect the Plug C connector to Position 2.</div> <div></div>
13	Install the top cover of 3000 Micro GC.
14	<div>Turn the 3000 Micro GC power switch, located on its front, to the on position. The switch will illuminate green to indicate that 3000 Micro GC is on.</div> <div>Wait three minutes for the instrument to load.</div>

This page is intentionally blank.